

**NASA/TM-2003-**

**Ocean Optics Protocols For Satellite Ocean Color Sensor  
Validation, Revision 5, Volume V:**

**Biogeochemical and Bio-Optical Measurements and Data Analysis  
Protocols**

*James L. Mueller, Giuletta S. Fargion and Charles R. McClain, Editors  
J. L. Mueller, R. R. Bidigare, C. Trees, W. M. Balch, J. Dore, D.T. Drapeau, D. Karl, L. Van  
Heukelem, and J. Perl Authors.*

National Aeronautical and  
Space administration

**Goddard Space Flight Space Center**  
Greenbelt, Maryland 20771

December 2003

**NASA/TM-2003-**

**Ocean Optics Protocols For Satellite Ocean Color Sensor  
Validation, Revision 5, Volume V:**

**Biogeochemical and Bio-Optical Measurements and Data  
Analysis Protocols**

*James L. Mueller, CHORS, San Diego State University, San Diego, California  
Giulietta S. Fargion, Science Applications International Corporation, Beltsville, Maryland  
Charles R. McClain, Goddard Space Flight Center, Greenbelt, Maryland*

*C. C. Trees and J. Perl, CHORS, San Diego State University, San Diego, California  
R. R. Bidigare, D. M. Karl and J. Dore, Dept. of Oceanography, University of Hawaii, Hawaii  
W. M. Balch and D. T. Drapeau, Bigelow Laboratory for Ocean Sciences, E. Boothbay Harbor, Maine  
L. Van Heukelem, University of Maryland Center for Environmental Science, Maryland*

National Aeronautical and  
Space administration

**Goddard Space Flight Space Center**  
Greenbelt, Maryland 20771

December 2003

## *Preface To Revision 5*

This document stipulates protocols for measuring bio-optical and radiometric data for the Sensor Intercomparison and Merger for Biological and Interdisciplinary Oceanic Studies (SIMBIOS) Project activities and algorithm development. The document is organized into 6 separate volumes, and in Revision 5, Volume VI is divided into 2 parts. Revision 5 consists of a new version of Volume V (Biogeochemical and Bio-Optical Properties) that supercedes and replaces Volume V (Revision 4), and new additions to Volume VI (Special Topics) are issued as Part 2 of that volume. The currently effective ocean optics protocol volumes, as of Revision 5, are:

### **Ocean Optics Protocols for Satellite Ocean Color Sensor Validation**

**Volume I: Introduction, Background and Conventions (Rev. 4)**

**Volume II: Instrument Specifications, Characterization and Calibration (Rev. 4)**

**Volume III: Radiometric Measurements and Data Analysis Methods (Rev. 4)**

**Volume IV: Inherent Optical Properties: Instruments, Characterization, Field Measurements and Data Analysis Protocols (Rev. 4 and Erratum 1 dated 28 Aug. 2003)**

**Volume V: Biogeochemical and Bio-Optical Measurements and Data Analysis Methods (Rev. 5)**

**Volume VI: Special Topics in Ocean Optics Protocols and Appendices (Rev. 4)**

**Volume VI, Part 2: Special Topics in Ocean Optics Protocols, Part 2 (Rev. 5)**

**Volume V (Revision 5):** This volume is issued as a complete replacement for **Volume V (Revision 4)**. The overview chapter (Chapter 1) briefly reviews biogeochemical and bio-optical measurements, and points to literature covering methods for measuring these variables. Detailed protocols for HPLC measurement of phytoplankton pigment concentrations are given in Chapter 2, and the Revision 5 version incorporates the **Erratum** issued in June 2003 to modify the HPLC protocols related to water retention by GF/F filters. Chapter 3 gives protocols for Fluorometric measurement of chlorophyll *a* concentration, and is carried over unchanged from Revision 4. Chapter 4 is a new addition which describes protocols for determining backscattering by Coccolithophorids and detached Coccoliths.

**Volume VI, Part 2 (Revision 5):** This volume supplements the 5 chapters of Volume VI (Rev. 4), adding two new “Special Topics” chapters:

- Chapter 6 briefly reviews recent progress in protocols for instrument self shading corrections to in-water upwelled radiance measurements;
- Chapter 7 reviews recent advances in radiometric characterization and measurement methods that are directly relevant to ocean color remote sensing and validation of satellite ocean color sensors.

This technical report is not meant as a substitute for scientific literature. Instead, it will provide a ready and responsive vehicle for the multitude of technical reports issued by an operational Project. The contributions are published as submitted, after only minor editing to correct obvious grammatical or clerical errors.

## *Preface To Revision 4*

This document stipulates protocols for measuring bio-optical and radiometric data for the Sensor Intercomparison and Merger for Biological and Interdisciplinary Oceanic Studies (SIMBIOS) Project activities and algorithm development. The document is organized into 6 separate volumes as:

### **Ocean Optics Protocols for Satellite Ocean Color Sensor Validation, Revision 4**

**Volume I: Introduction, Background and Conventions**

**Volume II: Instrument Specifications, Characterization and Calibration**

**Volume III: Radiometric Measurements and Data Analysis Methods**

**Volume IV: Inherent Optical Properties: Instruments, Characterization, Field Measurements and Data Analysis Protocols**

**Volume V: Biogeochemical and Bio-Optical Measurements and Data Analysis Methods**

**Volume VI: Special Topics in Ocean Optics Protocols and Appendices**

The earlier version of *Ocean Optics Protocols for Satellite Ocean Color Sensor Validation, Revision 3* (Mueller and Fargion 2002, Volumes 1 and 2) is entirely superseded by the seven Volumes of Revision 4 listed above.

The new multi-volume format for publishing the ocean optics protocols is intended to allow timely future revisions to be made reflecting important evolution of instruments and methods in some areas, without reissuing the entire document. Over the years, as existing protocols were revised, or expanded for clarification, and new protocol topics were added, the ocean optics protocol document has grown from 45pp (Mueller and Austin 1992) to 308pp in Revision 3 (Mueller and Fargion 2002). This rate of growth continues in Revision 4. The writing and editorial tasks needed to publish each revised version of the protocol manual as a single document has become progressively more difficult as its size increases. Chapters that change but little, must nevertheless be rewritten for each revision to reflect relatively minor changes in, *e.g.*, cross-referencing and to maintain self-contained consistency in the protocol manual. More critically, as it grows bigger, the book becomes more difficult to use by its intended audience. A massive new protocol manual is difficult for a reader to peruse thoroughly enough to stay current with and apply important new material and revisions it may contain. Many people simply find it too time consuming to keep up with changing protocols presented in this format - which may explain why some relatively recent technical reports and journal articles cite Mueller and Austin (1995), rather than the then current, more correct protocol document. It is hoped that the new format will improve community access to current protocols by stabilizing those volumes and chapters that do not change significantly over periods of several years, and introducing most new major revisions as new chapters to be added to an existing volume without revision of its previous contents.

The relationships between the Revision 4 chapters of each protocol volume and those of Revision 3 (Mueller and Fargion 2002), and the topics new chapters, are briefly summarized below:

**Volume I:** This volume covers perspectives on ocean color research and validation (Chapter 1), fundamental definitions, terminology, relationships and conventions used throughout the protocol document (Chapter 2), requirements for specific *in situ* observations (Chapter 3), and general protocols for field measurements, metadata, logbooks, sampling strategies, and data archival (Chapter 4). Chapters 1, 2 and 3 of Volume I correspond directly to Chapters 1, 2 and 3 of Revision 3 with no substantive changes. Two new variables, Particulate Organic Carbon (POC) and Particle Size Distribution (PSD) have been added to Tables 3.1 and 3.2 and the related discussion in Section 3.4; protocols covering these measurements will be added in a subsequent revision to Volume V (see below). Chapter 4 of Volume I combines material from Chapter 9 of Revision 3 with a brief summary of SeaBASS policy and archival requirements (detailed SeaBASS information in Chapter 18 and Appendix B of Revision 3 has been separated from the optics protocols).

**Volume II:** The chapters of this volume review instrument performance characteristics required for *in situ* observations to support validation (Chapter 1), detailed instrument specifications and underlying rationale (Chapter 2) and protocols for instrument calibration and characterization standards and methods (Chapters 3 through 5). Chapters 1 through 5 of Volume II correspond directly to Revision 3 chapters 4 through 8, respectively, with only minor modifications.

**Volume III:** The chapters of this volume briefly review methods used in the field to make the *in situ* radiometric measurements for ocean color validation, together with methods of analyzing the data (Chapter 1), detailed measurement and data analysis protocols for in-water radiometric profiles (Chapter 2), above water measurements of remote sensing reflectance (Chapter III-3), determinations of exact normalized water-leaving

radiance (Chapter 4), and atmospheric radiometric measurements to determine aerosol optical thickness and sky radiance distributions (Chapter 5). Chapter 1 is adapted from relevant portions of Chapter 9 in Revision 3. Chapter 2 of Volume III corresponds to Chapter 10 of Revision 3, and Chapters 3 through 5 to Revision 3 Chapters 12 through 14, respectively. Aside from reorganization, there are no changes in the protocols presented in this volume.

**Volume IV:** This volume includes a chapter reviewing the scope of inherent optical properties (IOP) measurements (Chapter 1), followed by 4 chapters giving detailed calibration, measurement and analysis protocols for the beam attenuation coefficient (Chapter 2), the volume absorption coefficient measured *in situ* (Chapter 3), laboratory measurements of the volume absorption coefficients from discrete filtered seawater samples (Chapter 4), and *in situ* measurements of the volume scattering function, including determinations of the backscattering coefficient (Chapter 5). Chapter 4 of Volume IV is a slightly revised version of Chapter 15 in Revision 3, while the remaining chapters of this volume are entirely new contributions to the ocean optics protocols. These new chapters may be significantly revised in the future, given the rapidly developing state-of-the-art in IOP measurement instruments and methods.

**Volume V:** The overview chapter (Chapter 1) briefly reviews biogeochemical and bio-optical measurements, and points to literature covering methods for measuring these variables; some of the material in this overview is drawn from Chapter 9 of Revision 3. Detailed protocols for HPLC measurement of phytoplankton pigment concentrations are given in Chapter 2, which differs from Chapter 16 of Revision 3 only by its specification of a new solvent program. Chapter 3 gives protocols for Fluorometric measurement of chlorophyll *a* concentration, and is not significantly changed from Chapter 17 of Revision 3. New chapters covering protocols for measuring, Phycoerythrin concentrations, Particle Size Distribution (PSD) and Particulate Organic Carbon (POC) concentrations are likely future additions to this volume.

**Volume VI:** This volume gathers chapters covering more specialized topics in the ocean optics protocols. Chapter 1 introduces these special topics in the context of the overall protocols. Chapter 2 is a reformatted, but otherwise unchanged, version of Chapter 11 in Revision 3 describing specialized protocols used for radiometric measurements associated with the Marine Optical Buoy (MOBY) ocean color vicarious calibration observatory. The remaining chapters are new in Revision 4 and cover protocols for radiometric and bio-optical measurements from moored and drifting buoys (Chapter 3), ocean color measurements from aircraft (Chapter 4), and methods and results using LASER sources for stray-light characterization and correction of the MOBY spectrographs (Chapter 5). In the next few years, it is likely that most new additions to the protocols will appear as chapters added to this volume. This volume also collects appendices of useful information. Appendix A is an updated version of Appendix A in Revision 3 summarizing characteristics of past, present and future satellite ocean color missions. Appendix B is the List of Acronyms used in the report and is an updated version of Appendix C in Revision 3. Similarly, Appendix C, the list of Frequently Used Symbols, is an updated version of Appendix D from Rev. 3. The SeaBASS file format information given in Appendix B of Revision 3 has been removed from the protocols and is promulgated separately by the SIMBIOS Project.

In the Revision 4 multi-volume format of the ocean optics protocols, Volumes I, II and III are unlikely to require significant changes for several years. The chapters of Volume IV may require near term revisions to reflect the rapidly evolving state-of-the-art in measurements of inherent optical properties, particularly concerning instruments and methods for measuring the Volume Scattering Function of seawater. It is anticipated that new chapters will be also be added to Volumes V and VI in Revision 5 (2003).

This technical report is not meant as a substitute for scientific literature. Instead, it will provide a ready and responsive vehicle for the multitude of technical reports issued by an operational Project. The contributions are published as submitted, after only minor editing to correct obvious grammatical or clerical errors.

*Table of Contents*

<b>CHAPTER 1.....</b>	<b>1</b>
<i>OVERVIEW OF BIOGEOCHEMICAL MEASUREMENTS AND DATA ANALYSIS IN OCEAN COLOR RESEARCH</i>	
1.1 INTRODUCTION .....	1
1.2 PHYTOPLANKTON PIGMENT CONCENTRATIONS .....	1
<i>High Performance Liquid Chromatography (HPLC) Measurements and Analysis</i> .....	1
<i>Fluorometric Measurement of Chlorophyll a Concentration</i> .....	1
<i>Phycocerythrin and other Phycobiliproteins</i> .....	2
1.3 IN SITU CHLOROPHYLL A FLUORESCENCE .....	2
1.4 SUSPENDED PARTICLES .....	3
<i>Suspended Particulate Matter</i> .....	3
<i>Particulate Organic Carbon and Particulate Organic Nitrogen</i> .....	3
<i>Particle Size Distributions</i> .....	3
<i>Coccolith Concentrations</i> .....	3
1.5 FUTURE DIRECTIONS .....	3
REFERENCES .....	4
<b>CHAPTER 2.....</b>	<b>5</b>
<i>HPLC PHYTOPLANKTON PIGMENTS: SAMPLING, LABORATORY METHODS, AND QUALITY ASSURANCE PROCEDURES</i>	
2.1 INTRODUCTION .....	5
2.2 SAMPLING PROTOCOLS FOR PHYTOPLANKTON PIGMENTS .....	6
<i>Water Samples</i> .....	6
<i>Filtration</i> .....	6
<i>Sample Handling and Storage</i> .....	7
<i>Recordkeeping</i> .....	8
2.3 LABORATORY METHODS FOR HPLC PHYTOPLANKTON PIGMENT ANALYSIS .....	8
<i>Internal Standard and Solvent Preparation</i> .....	8
<i>Extraction</i> .....	8
<i>Apparatus</i> .....	9
<i>HPLC Eluants and Gradient Programs</i> .....	9
<i>Determination of Algal Chlorophyll and Carotenoid Pigments by HPLC (Wright et al. 1991):</i> .....	9
2.4 QUALITY ASSURANCE PROCEDURES .....	11
2.5 PROTOCOL STATUS AND FUTURE DIRECTIONS FOR RESEARCH .....	12
REFERENCES .....	12
<b>CHAPTER 3.....</b>	<b>15</b>
<i>FLUOROMETRIC CHLOROPHYLL A: SAMPLING, LABORATORY METHODS, AND DATA ANALYSIS PROTOCOLS</i>	
3.1 INTRODUCTION .....	15
3.2 SAMPLE ACQUISITION AND STORAGE .....	16
<i>Filtration</i> .....	16
<i>Sample Handling, and Storage</i> .....	17
<i>Recordkeeping</i> .....	17
3.3 LABORATORY METHODS FOR FLUOROMETRIC DETERMINATION OF CHL. A AND PHEOPIGMENT CONCENTRATIONS .....	17
<i>Fluorometer Calibrations</i> .....	18
<i>Solvent Preparation</i> .....	19
<i>Extraction</i> .....	19
<i>Measurement</i> .....	19
3.4 IN SITU CHLOROPHYLL A FLUORESCENCE PROFILES .....	20
3.5 PROTOCOL STATUS AND FUTURE DIRECTIONS FOR RESEARCH .....	20
REFERENCES .....	24

**CHAPTER 4..... 27**

*BACKSCATTERING BY COCCOLITHOPHORIDS AND COCCOLITHS: SAMPLE PREPARATION, MEASUREMENT AND ANALYSIS PROTOCOLS*

4.1 INTRODUCTION .....	27
4.2 VOLUME SCATTERING FUNCTION INSTRUMENTS AND CALIBRATION .....	28
4.3 DISCRETE WATER SAMPLES: SAMPLE ACQUISITION, TREATMENT, AND VSF MEASUREMENTS.....	31
<i>Sample Acquisition</i> .....	31
<i>VSF Measurements of Untreated Seawater Samples</i> .....	31
<i>Sample Acidification to Dissolve Particulate CaCO<sub>3</sub></i> .....	31
<i>VSF Measurements of Acidified Samples</i> .....	32
<i>Quality Control and Uncertainty</i> .....	32
4.4 CONTINUOUS FLOW-THROUGH MEASUREMENTS ALONG SHIP TRACKS: SAMPLE ACQUISITION, TREATMENT, AND VSF MEASUREMENTS .....	32
<i>Flow-Through Sampling Apparatus and Debubbler</i> .....	32
<i>In Line Acidification and Flushing Method</i> .....	32
<i>Measurement Cycles: Untreated Sample VSF – Acidified Sample VSF – System Flushing</i> .....	32
<i>Quality Control</i> .....	33
4.5 ANCILLARY MEASUREMENTS.....	33
<i>Coccolith and Coccolithophore Enumeration</i> .....	33
<i>Particulate Inorganic Carbon</i> .....	33
4.6 DATA ANALYSIS METHODS .....	34
4.7 DISCUSSION: PROTOCOL STATUS AND FUTURE DIRECTIONS .....	34
ACKNOWLEDGEMENTS.....	35
REFERENCES .....	35

## Chapter 1

# Overview of Biogeochemical Measurements and Data Analysis in Ocean Color Research

James L. Mueller

*Center for Hydro-Optics and Remote Sensing, San Diego State University, California*

### 1.1 INTRODUCTION

A total of 9 biogeochemical and bio-optical observations are listed in Tables 3.1 and 3.2 (Vol. I). Phytoplankton pigment concentrations determined by the HPLC method, and fluorometric chlorophyll *a* and pheopigment concentrations are **required** measurements for which detailed protocols are described in Vol. I, Chapters 2 and 3, respectively. Observation of chlorophyll *a* fluorescence intensity *in situ* is listed as highly desired, and protocols for its measurement and data analysis are also included in Chapter 3. Six additional biogeochemical observations are listed as **specialized measurements**. These include concentrations of Phycobiliprotein (Phycocerythrin), and suspended particulate measurements including *Coccolith* concentrations, total Suspended Particulate Matter (SPM), Particulate Organic Carbon (POC), Particulate Organic Nitrogen (PON, and Particle Size Distribution (PSD). Methods for measurement and data analysis for backscattering by *Coccolithophorids* and detached *Coccoliths* are described in Chapter 4, a new addition in Revision 5. Methods of measurement and data analysis for these specialized observations, most of which are related to applications of ocean color image data to ocean process studies, are reviewed briefly in the present chapter.

### 1.2 PHYTOPLANKTON PIGMENT CONCENTRATIONS

#### *High Performance Liquid Chromatography (HPLC) Measurements and Analysis (Chapter 2)*

Mueller and Austin (1995) simply adopted the JGOFS HPLC protocols for measuring phytoplankton pigment concentrations by reference (UNESCO 1994), and supplemented them with some brief instructions on sampling and sample handling procedures. Although this approach embraced protocol documentation describing a complete methodology, and represented a community consensus, the lack of a comprehensive end-to-end protocol statement has proved to be a source of confusion and debate within the ocean color community. Furthermore, the JGOFS protocols (UNESCO 1994) specified that pigment concentrations should be reported in units of pigment mass per mass of seawater ( $\text{ng Kg}^{-1}$ ), rather than in units of pigment mass per volume of seawater (either  $\mu\text{g L}^{-1}$ , or  $\text{mg m}^{-3}$ ). The use of volumetric concentrations is critical because radiative transfer in the ocean, and absorption by pigments, are volumetric processes. One could use the mass concentration values preferred by JGOFS, but it would be essential to supplement them with densities computed from CTD data, and make the conversion to volumetric concentrations. Therefore, a complete set of protocols for HPLC measurement of phytoplankton pigment concentrations was added as Chapter 13 of Revision 2.0 to the Ocean Optics Protocols (Fargion and Mueller 2000), updated as Chapter 16 of Revision 3 (Mueller and Fargion 2002), and updated here again as Chapter 2 of the present volume. Chapter 2 provides complete protocols for obtaining water samples, filtering them, freezing the filtered samples in liquid nitrogen, sample handling and storage, extraction, HPLC calibrations and measurements, data analysis and quality control. A new HPLC solvent program in Chapter 2 replaces that specified in the previous version of the protocols (Bidegare *et al.* 2002).

#### *Fluorometric Measurement of Chlorophyll *a* Concentration (Chapter 3)*

For reasons similar to those described above for HPLC pigment measurements, it was decided that the protocols for fluorometric measurement of the concentrations of chlorophyll *a* and phaeopigments were too briefly abstracted in Mueller and Austin (1995). Therefore, new detailed protocols for this measurement were added as Chapter 14 to Revision 2 (Fargion and Mueller 2000), updated as Chapter 17 of Revision 3 (Mueller and Fargion 2002), and reproduced here without significant change as Chapter 3. Chapter 3 provides complete protocols for obtaining water

samples, filtering them, freezing the filtered samples in liquid nitrogen, sample handling and storage, extraction, fluorometer calibrations and measurements, data analysis and quality control.

In addition, Chapter 3 discusses geographic and temporal variability in the relationship between fluorometric chlorophyll concentrations and combined concentrations of total chlorophyll pigments determined by the HPLC methods (Chapter 2). It is both easier and less expensive to measure chlorophyll *a* and pheopigment concentrations using the fluorometric method, which has the added advantage of allowing shipboard analyses at sea during lengthy cruises. When these data are used for remote sensing algorithm development or validation, however, regional and temporal (*i.e.* cruise-to-cruise) dispersions and/or biases may be introduced unless the fluorometric data are first statistically adjusted (on a local basis) to agree with HPLC determinations of the concentration of total chlorophylls. A cost-effective strategy is to acquire, on each cruise, a majority of filtered pigment samples for fluorometric chlorophyll *a* and pheopigment analysis, supplemented by a smaller number of replicate samples for HPLC pigment analysis. The HPLC replicates should provide a representative distribution over geographic location, depth and time during a cruise, and will be used to determine a local regression relationship between the two measurements. This approach is now required for pigment data submitted for SeaBASS archival and SIMBIOS validation analysis.

#### *Phycocerythrin and other Phycobiliproteins*

$R_{RS}(\lambda)$  may be enhanced by fluorescence by phycocerythrin (PE) in a band near 565 nm (*e.g.* Hoge *et al.* 1998; Wood *et al.* 1999). The detection from aircraft of laser-induced phycocerythrin fluorescence is already well established (Hoge *et al.* 1998). It is more difficult to detect and quantify solar induced phycocerythrin fluorescence, but some work has been done in that area as well (Morel *et al.* 1993; Morel 1997; Hoge *et al.* 1999; Subramaniam *et al.* 1999).

Various phycocerythrins differ from one another in chromophore composition. All phycocerythrins contain phycocerythrobilin chromophores [PEB; maximum  $a(\lambda)$  near  $\lambda \sim 550$  nm]; many others also contain phycocourobilin chromophores [PUB; maximum  $a(\lambda)$  near  $\lambda \sim 500$  nm], which extends the range of wavelengths absorbed by the pigment molecule into the blue regions of the spectrum. The ratio of PUB:PEB chromophores in the PE pigments synthesized by different *Synechococcus* strains greatly affects the absorption spectrum of the whole cells (Wood *et al.* 1985). Clearly, the dependence of  $a(\lambda)$  on the PUB:PEB ratio of phycocerythrin will affect also  $R_{RS}(\lambda)$  in water masses dominated by cyanobacteria. The PUB:PEB ratio for the PE in a given water mass may be characterized using scanning fluorescence spectroscopy (Wood *et al.*, 1999; Wyman, 1992).

The measurement of phycocerythrin is not yet as routine, nor as accurate, as the measurements of chlorophylls and carotenoids. The techniques introduced by Stewart and Farmer (1984) work well for measuring biliproteins in freshwater and estuarine species, but are less successful for natural populations of marine species. Wyman (1992) reported a linear relationship between the *in vivo* fluorescence emission intensity of PE measured in the presence of glycerol and the PE content of *Synechococcus* strain WH7803. Scanning spectral fluorescence measurements have been used to estimate PE concentration of extracted bulk samples (Vernet *et al.*, 1990). Nevertheless, there are few direct measurements of separated PE proteins from natural samples. High Performance Capillary Electrophoresis (HPCE) is a powerful analytical tool currently used in clinical, biochemical, pharmaceutical, forensic, and environmental research. In HPCE, high voltages (typically 10-30 KV) are used to separate molecules rapidly in narrow-bore (25-100  $\mu$ m), fused-silica capillaries based on differences in the charge-to-mass ratio of the analytes. HPCE is an automated analytical separation system with reduced analysis times and on-line quantification of compounds, ideally suited to the separation and quantification of water-soluble proteins (like phycobilins) from seawater. HPCE methods for separation analyses of phycocerythrin from cyanobacterial cultures and natural samples are currently under development and may be included in a future revision to the ocean optics protocols (C. Kinkade, Pers. Comm.).

### **1.3 IN SITU CHLOROPHYLL *a* FLUORESCENCE**

Protocols for measuring and analyzing profiles of *in situ* fluorescence by chlorophyll *a*,  $F(z)$  (Table 3.1 in Chapter 3, Volume I) are described in Chapter 3. When measured together with  $c(z,660)$  profiles (Chapter 2, Volume IV), the structure of  $F(z)$  provides valuable guidance for selecting depths of water samples, analyses of structure in  $K(z,\lambda)$  derived from radiometric profiles, and various aspects of quality control analysis. It is often useful to digitally record one-minute averages of  $F(z, lat, lon)$  in water pumped from a near-surface depth ( $z \sim 3$  m) to measure horizontal variability while underway steaming between stations, especially in water masses where

mesoscale and sub-mesoscale variability is strong (Section 4.2, Chapter 4, Vol. I). If supplemented by frequent fluorometric chlorophyll *a* samples filtered from the flow-through system, the alongtrack profile of  $F(z, lat, lon)$  can be “calibrated” in units of chlorophyll *a* concentration ( $\text{mg m}^{-3}$ ).

## 1.4 SUSPENDED PARTICLES

### *Suspended Particulate Matter*

All total suspended particulate material (SPM) dry weight ( $\text{mg L}^{-1}$ ) will be determined gravimetrically as outlined in Strickland and Parsons (1972)<sup>1</sup>. In general, samples are filtered through preweighed  $0.4 \mu\text{m}$  polycarbonate filters. The filters are washed with three 2.5 mL - 5.0 mL aliquots of DIW and immediately dried, either in an oven at  $75^\circ \text{C}$ , or in a dessicator. The filters are then reweighed in a laboratory, back on shore, using an electrobalance with at least seven digits of precision.

### *Particulate Organic Carbon and Particulate Organic Nitrogen*

Protocols for measuring concentrations in seawater of Particulate Organic Carbon (POC) and Particulate Organic Nitrogen (PON), as specified for JGOFS (UNESCO 1994, Chapter 15), are also adopted here. The units of POC and PON are  $\mu\text{g C Kg}^{-1}$  and  $\mu\text{g N Kg}^{-1}$ , respectively. Therefore, it is mandatory that each of these measurements be accompanied by Conductivity, Temperature and Pressure measurements so that the density of seawater [ $\text{Kg m}^{-3}$ ] may be calculated.

### *Particle Size Distributions*

Particle size distributions can potentially provide important information about the shape of the volume scattering function, which strongly influences the bidirectional aspects of remote-sensing reflectance (Chapter 4 of Volume III and, *e.g.*, Morel and Gentili 1996). Particle size distributions have been measured for many years using Coulter Counters and related to IOP, including  $c(\lambda)$  (*e.g.* Kitchen *et al.* 1982). More recently, several investigators have used the Spectrix Particle Size Analyzer to measure particle size distributions (see, *e.g.*, Chapter 2, Vol. VI). Protocols for measurements and analyses of particle size distributions are not included in this version of the ocean optics protocols, but should be written and added to a future revision of this protocol volume.

### *Coccolith Concentrations (Chapter 4)*

Concentrations of coccoliths, calcium carbonate ( $\text{CaCO}_3$ ) platelets detached from coccolithophorids (*sp.*), are measured as cell counts [number density per unit volume] using a microscope with polarization optics (Balch *et al.* 1991). An epifluorescence microscope is used to count plated and naked intact cells, before and after the coccoliths are dissolved by acidification. Also measured, before and after acidification, are the Volume Scattering Function (VSF) values at many angles, from which the volume specific backscattering coefficient for coccoliths is determined by subtraction (Voss *et al.* 1998). These methods are described in Chapter 4, which is a new addition to the protocols in Revision 5.

## 1.5 FUTURE DIRECTIONS

Future additions to this volume may include chapters providing detailed protocols for Phycoerythrin measurement and data analysis, and methods for measurement and analyses of SPM, PSD, and the organic suspended particulate fractions POC and PON.

---

<sup>1</sup> In some previous versions of the Ocean Optics Protocols (Mueller and Austin 1992, 1995; Fargion and Mueller 2000), it was incorrectly stated that suitable protocols were part of the JGOFS core measurements protocols (UNESCO 1994). The JGOFS protocols do not include SPM measurements of the type specified here.

## REFERENCES

- Balch, W.M., P.M. Holligan, S.G. Ackleson, and K.J. Voss, 1991: Biological and optical properties of mesoscale coccolithophore blooms in the Gulf of Maine. *Limnol. Oceanogr.*, **36**: 629-643.
- Bidegare, R.R., L. Van Heukelem, and C.C. Trees, 2002: HPLC phytoplankton pigments: sampling, laboratory methods, and quality assurance procedures. Chapter 17 in *Ocean Optics Protocols for Satellite Ocean Color Sensor Validation, Revision 3*, Mueller J.L. and G.S. Fargion [Eds.], NASA TM 2002-210004, NASA Goddard Space Flight Center, Greenbelt, Maryland, pp258-268.
- Fargion, G.S. and J.L. Mueller, 2000: *Ocean Optics Protocols for Satellite Ocean Color Sensor Validation, Revision 2*, NASA TM 2001-209955, NASA Goddard Space Flight Center, Greenbelt, Maryland, 184 pp.
- Hoge, F. E., C. W. Wright, P. E. Lyon, R. N. Swift, and J. Yungel, 1999: Satellite retrieval of the absorption coefficient of phytoplankton phycoerythrin pigment: Theory and feasibility status. MODIS ATBD doc. 27.
- Kitchen, J.C., J.R.V. Zaneveld and H. Pak, 1982: Effect of particle size distribution and chloropyll content on beam attenuation spectra. *Appl. Opt.*, **21**: 3913-3918.
- Morel, A. 1997: Consequences of a *Synechococcus* bloom upon the optical properties of oceanic (case 1) waters. *Limnol. Oceanogr.*, **42**: 1746-1754.
- Morel, A., Y.H. Ahn, F. Partensky, D. Vaultot, and H. Claustre. 1993: *Prochlorococcus* and *Synechococcus*: A comparative study of their optical properties in relation to their size and pigmentation. *J. Mar. Res.*, **51**: 617-647.
- Mueller, J.L., and R.W. Austin, 1995: Ocean Optics Protocols for SeaWiFS Validation, Revision 1. *NASA Tech. Memo. 104566, Vol. 25*, S.B. Hooker, E.R. Firestone and J.G. Acker, Eds., NASA Goddard Space Flight Center, Greenbelt, Maryland, 67 pp.
- Mueller, J.L. and G.S. Fargion, 2002: *Ocean Optics Protocols for Satellite Ocean Color Sensor Validation, Revision 3*, NASA TM 2002-210004, NASA Goddard Space Flight Center, Greenbelt, Maryland, 184 pp.
- Stewart, D.E. and F.H. Farmer. 1984: Extraction, identification, and quantification of phycobiliprotein pigments from phototrophic plankton. *Limnol. Oceanogr.*, **29**: 392-397.
- Strickland, J.D.H., and T.R. Parsons, 1972: *A Practical Handbook of Sea Water Analysis*, Fisheries Research Board of Canada, 310 pp.
- Subramaniam, A., E. J. Carpenter, and P. G. Falkowski, 1999: Bio-optical properties of the marine diazotrophic cyanobacteria *Trichodesmium* spp. II. A reflectance model for remote sensing. *Limnol. Oceanogr.*, **44**: 618-627.
- UNESCO, 1994: Protocols for the Joint Global Ocean Flux Study (JGOFS) Core Measurements, Manuals and Guides **29**: 170pp
- Vernet, M., B.G. Mitchell, and O. Holm-Hansen: 1990: Adaptation of *Synechococcus in situ* determined by variability in intracellular phycoerythrin-543 at a coastal station off the Southern California coast, USA. *Mar. Ecol. Prog. Ser.*, **63**: 9-16.
- Voss, K.J., W.M. Balch and K.A. Kilpatrick, 1998,. Scattering and attenuation properties of *Emiliania huxleyi* cells and their detached coccoliths. . *Limnol. Oceanogr.*, **43**(5): 870-876.
- Wood, A.M., P.K. Horan, K. Muirhead, D. Phinney, C.M. Yentsch, and J.M. Waterbury, 1985: Discrimination between types of pigments in marine *Synechococcus* by scanning spectroscopy, epifluorescence microscopy, and flow cytometry. *Limnol. and Oceanogr.*, **30**: 1303-1315.
- Wood, A.M., M. Lipsen and P. Coble, 1999: Fluorescence based characterization of phycoerythrin-containing cyanobacterial communities in the Arabian Sea during the Notheast and early Southwest Monsoon (1994-1995). *Deep-Sea Res. II*, **46**: 1769-1790.
- Wyman, M. 1992: An in vivo method for the estimation of phycoerythrin concentrations in marine cyanobacteria (*Synechococcus* spp.). *Limnol. Oceanogr.*, **37**: 1300-1306.

## Chapter 2

# HPLC Phytoplankton Pigments: Sampling, Laboratory Methods, and Quality Assurance Procedures

Robert R. Bidigare<sup>1</sup>, Laurie Van Heukelem<sup>2</sup> Charles C. Trees<sup>3</sup> and Jason Perl<sup>3</sup>

<sup>1</sup> *Department of Oceanography, University of Hawaii, Hawaii*

<sup>2</sup> *Horn Point Environmental Laboratory, University of Maryland, Maryland*

<sup>3</sup> *Center for Hydro-Optics and Remote Sensing, San Diego State University, California*

### 2.1 INTRODUCTION

Marine phytoplankton utilize chlorophyll *a* as their major light harvesting pigment for photosynthesis. Other accessory pigment compounds, such as chlorophylls *b* and *c*, carotenoids and phycobiliproteins, also play a significant role in photosynthesis by extending the organism's optical collection window, thereby improving absorption efficiencies and adaptation capabilities. The important chlorophyll degradation products found in the aquatic environment are the chlorophyllides, phaeophorbides, and phaeophytins. The presence, or absence, of the various photosynthetic pigments is used to separate the major algal groups, and to map the chemotaxonomic composition of phytoplankton in the oceans.

The unique optical properties of chlorophyll *a* have been used to develop spectrophotometric (Jeffrey and Humphrey 1975) and fluorometric (Holm-Hansen *et al.* 1965) measurement techniques. With the commercial availability of fluorometers for routine measurements of chlorophyll *a*, this pigment became a universal parameter in biological oceanography for estimating phytoplankton biomass and productivity. These optical methods can significantly under- or overestimate chlorophyll *a* concentrations, because of the overlap of the absorption and fluorescence bands of co-occurring chlorophylls *b* and *c*, chlorophyll degradation products, and accessory pigments (Trees *et al.* 1985; Smith *et al.* 1987; Hoepffner and Sathyendranath 1992; Bianchi *et al.* 1995; Tester *et al.* 1995).

The application of HPLC to phytoplankton pigment analysis has lowered the uncertainty for measuring chlorophyll *a* and pheopigments, as well as the accessory pigments, since compounds are physically separated and individually quantified. HPLC has provided oceanographers with a powerful tool for studying the processes affecting the phytoplankton pigment pool. Pigment distribution is useful for quantitative assessment of phytoplankton community composition, phytoplankton growth rate and zooplankton grazing activity.

For low uncertainty determinations of chlorophylls *a*, *b*, and *c*, chlorophyll degradation products, and carotenoid pigments, HPLC techniques are recommended. It should be noted, however, that the reverse-phase C<sub>18</sub> HPLC method recommended by the Scientific Committee on Oceanographic Research (SCOR) (Wright *et al.* 1991) is not capable of separating monovinyl chlorophyll *a* from divinyl chlorophyll *a*, nor monovinyl chlorophyll *b* from divinyl chlorophyll *b*. This method, therefore, only provides concentration *estimates* for these co-eluting pigment pairs; methods for optically resolving monovinyl chlorophyll *a* and divinyl chlorophyll *a* are given below.

Divinyl chlorophyll *a*, the major photosynthetic pigment found in *Prochlorococcus*, accounts for 10 % to 60 % of the total chlorophyll *a* in subtropical and tropical oceanic waters (Goericke and Repeta 1993; Letelier *et al.* 1993; Andersen *et al.* 1996; Bidigare and Ondrusek 1996; Gibb *et al.* 2000). Divinyl chlorophyll *a* is spectrally different from *normal* (monovinyl) chlorophyll *a* and its presence results in a significant overestimation of total chlorophyll *a* concentration as determined by the conventional HPLC methods (Goericke and Repeta 1993; Letelier *et al.* 1993; Latasa *et al.* 1996). To avoid these errors, it is recommended that monovinyl and divinyl chlorophyll *a* be spectrally resolved, or chromatographically separated, to obtain an unbiased determination of total chlorophyll *a* for ground-truthing satellite ocean color algorithms and imagery. Total chlorophyll *a*, TChl *a*, is the sum of divinyl chlorophyll *a*, monovinyl chlorophyll *a*, chlorophyllide *a*, and chlorophyll *a* epimers and allomers. These co-eluting chlorophyll species can be resolved spectrally following C<sub>18</sub> HPLC chromatography (Wright *et al.* 1991) and quantified using dichromatic equations at 436 nm and 450 nm (Latasa *et al.* 1996). Alternatively, these two chlorophyll species can be separated chromatographically and individually quantified using C<sub>8</sub> HPLC techniques (see below).

The protocols specified below for HPLC pigment analyses follow closely those prescribed in the *JGOFS Core Measurement Protocols* (UNESCO 1994). Both sets of protocols include:

1. Use of Whatman GF/F glass fiber filters, approximately 0.7  $\mu\text{m}$  pore size;
2. Extraction in aqueous acetone; and
3. Calibration with standards.

The present protocols differ from the JGOFS protocols in one critical respect. Absorption of light in seawater, or any other medium, is a volumetric process, even though the volume absorption coefficient may vary with the density of the medium. For ocean color and optical analyses, therefore, the concentrations in seawater of all phytoplankton pigments shall be expressed in units of mass per unit volume of seawater ( $\mu\text{g L}^{-1}$  or  $\text{mg m}^{-3}$ ). This differs from the JGOFS protocols, which specify that concentrations in seawater of all phytoplankton pigments should be expressed in  $\text{ng Kg}^{-1}$ .

In addition to HPLC analyses, it is recommended that the standard fluorometric methodology used for measuring chlorophylls and pheopigments (Holm-Hansen *et al.* 1965, Strickland and Parson 1972) also be applied to the same extracted pigment samples used for HPLC analysis. Protocols for fluorometric measurements of chlorophyll *a* and pheopigments are given here in Chapter 3 of the present volume. For a more in depth review of guidelines for measuring phytoplankton pigments in oceanography see Jeffrey *et al.* (1997)

## 2.2 SAMPLING PROTOCOLS FOR PHYTOPLANKTON PIGMENTS

### *Water Samples*

Water samples should be taken using, *e.g.*, Niskin bottles at the site of, and simultaneously with, the surface in-water upwelled radiance and reflectance measurements, and at depth increments sufficient to resolve variability within at least the top optical depth. The  $K(z,\lambda)$ , profiles over this layer will be used to compute optically weighted, near-surface pigment concentration for bio-optical algorithm development (Gordon and Clark 1980).

When possible, samples should be acquired at several depths distributed throughout the upper 200 m of the water column [or in turbid water, up to seven diffuse attenuation depths, *i.e.*  $\ln(E(z,\lambda)/E(z,\lambda)_0)=7$ ], to provide a basis for relating fluorescence signals to pigment mass concentrations.

Samples should be filtered as soon as possible after collection. If processing must be delayed for more than an hour, hold the samples on ice, or in a freezer at 4°C, and protect them from exposure to light. For delays longer than several hours, the samples should be stored in liquid nitrogen. Use opaque sample bottles, because even brief exposure to light during sampling and/or storage might alter pigment values.

### *Filtration*

Whatman GF/F glass fiber filters, with approximately 0.7  $\mu\text{m}$  pore size, are preferred for removing phytoplankton from water. The glass fibers assist in breaking the cells during grinding, accommodate larger sample volumes, and do not form precipitates after acidification. Twenty-five mm diameter GF/F glass fiber filters should be used with vacuum (7-8 inches of mercury) or positive pressure (1-2 psi). Positive pressure filtration is recommended, because it filters larger volumes of water at reduced filtration times. The only problem with vacuum filtration is that unobservable air leaks may occur around the filtration holder, and as a result the pressure gradient across the filter is much less than what is indicated on the vacuum gauge. When positive filtration is used, any leakage around the filter holder results in observable dripping water.

Inert membrane filters, such as polyester filters, may be used when size fraction filtration is required. When this is done, it is recommended to also filter a replicate sample through a GF/F to determine the total concentration. Summing the various size-fractionated concentrations may not produce an accurate estimate of the total, because of the potential for cell disruption during filtration.

There has been an ongoing discussion of filter types and retention efficiencies for natural samples. Phinney and Yentsch (1985) showed the inadequacy of GF/F filters for retaining chlorophyll *a* in oligotrophic waters, as did Dickson and Wheeler (1993) for samples from the North Pacific. In response to Dickson and Wheeler (1993),

Chavez *et al.* (1995) compared samples collected in the Pacific Ocean using GF/F and 0.2  $\mu\text{m}$  membrane filters with small filtered volumes (100 mL to 540 mL). Their results showed a very close agreement between the two filter types, with GF/F filters having only a slightly positive 5 % bias.

Filtration volume can directly affect the retention efficiency for GF/F filters. Particles can be retained by filters through a variety of ways, such as filter sieving, filter adsorption, electrostatic and van der Waals attractions (Brock, 1983). When water flows through the pores of a Nuclepore filter, streamlines are formed that can align small particles longitudinally, with the result that cell diameter becomes important with these filters. It is known, on the other hand, that Whatman GF/F filters can retain particles much smaller than their rated pore size. Generally, at small volumes (100 mL to 300 mL) filter adsorption, and electrostatic and van der Waals attractions are important, whereas at larger volumes (>2,000 mL) sieving dominates. This has been tested in oligotrophic waters off Hawaii in which small (<500 mL) and large volumes (> 2 L to 4 L) retained similar amounts of chlorophyll *a* on the two types of filters, whereas for intermediate sample volumes the GF/F filters showed lower concentrations. During several cruises off the Hawaiian Islands, differences in retention efficiencies were found for GF/F filters to be a function of sample volume; large sample volumes (2 L and 4 L) retained about 18 % more chlorophyll *a* than replicate 1 L samples.

Filtration volumes are usually limited by the concentration of particles present in each sample. For HPLC analysis it is important to filter as large a volume as possible, so as to accurately measure most of the major pigments. A qualitative check to determine whether a large enough volume has been filtered is to count the number of accessory pigments (chlorophylls *b*, *c*<sub>1</sub>, *c*<sub>2</sub>, *c*<sub>3</sub>, and carotenoids) quantified, excluding chlorophyll degradation products (Trees *et al.* 2000). Most algal groups (excluding phycobiliprotein-containing groups) contain at least *four* HPLC-measurable accessory pigments (see Jeffrey *et al.* 1997). Therefore, pigment samples that do not meet this minimum accessory pigment criterion may have detection limit problems related to low signal-to-noise ratios for the HPLC detectors and/or inadequate concentration techniques (*e.g.* low filtration volumes). It is generally recommended that the following volumes be filtered for HPLC pigment analyses: 3 L to 4 L for oligotrophic waters, 1 L to 2 L for mesotrophic waters, and 0.5 L to 1 L for eutrophic waters.

It is recommended to not pre-filter seawater samples to remove large zooplankton and particles, because this practice may exclude pigment-containing colonial and chain-forming phytoplankton, such as diatoms and *Trichodesmium* sp. Forceps may be used to remove large zooplankton from the GF/Fs following filtration.

#### *Sample Handling and Storage*

Samples should be filtered as quickly as possible after collection and stored immediately in liquid nitrogen. Liquid nitrogen is the best method for storing samples with minimum degradation for short, as well as, longer storage times (*e.g.* 1 year). Placing samples in liquid nitrogen also assists in pigment extraction by weakening the cell wall and membrane during this rapid temperature change. Ultra-cold freezers (-90 °C) can be used for storage, although they have not been tested for longer than 60 days (Jeffrey *et al.* 1997). Conventional deep freezers should not be used for storing samples more than 20 hours before transferring them to an ultra-cold freezer, or liquid nitrogen. Again, storage of samples in liquid nitrogen immediately after filtration is the preferred method.

Samples should be folded in half with the filtered halves facing in. This eliminates problems of rubbing particles off the filter during placement in sample containers and storage.

It is strongly recommended to use aluminum foil wrappings for sample containers. This simple, but effective, container is both inexpensive and easy to use. Cut small pieces of heavy-duty aluminum foil into approximately 4 cm squares. Fold each piece in half, and using a fine-point permanent marker, write a short sample identifier (*e.g.* first letter of the cruise and a sequential sample number) on the foil. Writing on the folded foil, prior to placement of the filter, both avoids puncturing the foil with the marking pen, and improves the legibility of the sample identifier. Place the folded filter in the aluminum foil. Fold the three open sides to form an envelope that is only slightly larger than the folded filter (~3 cm x 1.5 cm).

The use of foil containers minimizes the size requirement of the storage container. It is also acceptable to use either cryogenic tubes, or HistoPrep tissue capsules, but they occupy more storage volume per sample, and they are more expensive than aluminum foil. If fluorometric analysis is to be done soon after collection, it is still recommended to place the samples in liquid nitrogen to assist in pigment extraction, and on removal from the liquid nitrogen to place them immediately in chilled 90 % acetone.

*Recordkeeping*

Information regarding sample identification should be logged in a laboratory notebook with the analyst's initials. For each filter sample record the sample identifier (as written on the sample container), station number for the cruise, water volume filtered ( $V_{\text{FILT}}$ ) in mL, and depth of the water sample, together with the date, time, latitude, and longitude of the bottle cast during which the sample was acquired.

## 2.3 LABORATORY METHODS FOR HPLC PHYTOPLANKTON PIGMENT ANALYSIS

*Internal Standard and Solvent Preparation*

In addition to daily calibration of the HPLC system with external standards, an internal standard (*e.g.* canthaxanthin) should be used to determine the extraction volume. It is important to verify that the internal standard employed is not a *naturally* occurring analyte in the field samples to be analyzed by HPLC. Canthaxanthin is recommended as an internal standard because it has a restricted distribution in ocean waters, and it is readily available in high purity from commercial sources. For additional background on the use of internal standards see Snyder and Kirkland (1979). The internal standard should be added to the sample prior to extraction and used to correct for the addition of GF/F filter-retained seawater and sample volume changes during extraction. When new external and internal standards are prepared they should be verified against previous standards and a standard reference solution if available. An internal standard with an HPLC peak removed from those of all the pigments, canthaxanthin, is added at a fixed concentration to the HPLC-grade acetone solvent used to extract the pigments from the filtered samples. A sample of canthaxanthin spiked solvent is injected into the HPLC system and its peak area  $A_{\text{STD}}^{\text{Cantha}}$  is recorded to provide a baseline internal standard for monitoring the solvent concentration in each extracted sample.

*Extraction*

Filters are removed from the liquid nitrogen, briefly thawed (~1 min), and placed in glass centrifuge tubes for extraction in acetone. Three mL HPLC-grade acetone is added to each tube, followed by the addition of a fixed volume of internal standard (typically 50  $\mu\text{L}$  canthaxanthin in acetone). Alternatively, canthaxanthin spiked HPLC-grade acetone solvent may be prepared in advance, in a batch large enough for all samples, and 3.00 mL is added to each tube in a single step. Since GF/F filters retain a significant amount of seawater following filtration (ca. 0.22 mL per 25 mm filter), the final acetone concentration in the pigment extracts is ~ 94 % (acetone:water, by volume); by measuring the canthaxanthin peak area  $A_{\text{STD}}^{\text{Cantha}}$  for each sample, the ratio  $A_{\text{STD}}^{\text{Cantha}} / A_{\text{Sample}}^{\text{Cantha}}$  may be used to adjust for sample to sample variations in the extraction volume. If a canthaxanthin internal standard is used to correct for volume differences caused by dilution, or evaporation, the volume extracted,  $V_{\text{extracted}}$ , is recorded as the unadjusted volume of solvent added during extraction. If an internal standard is not used, then  $V_{\text{extracted}}$  must be corrected for the water retained by the filter. The correction for seawater retention by a GF/F filter is approximately 0.22 mL, a value determined originally by Trees (1978) and recently confirmed in a laboratory experiment at CHORS by J. Perl.

Samples are disrupted by sonication, placed in a freezer, and allowed to extract at 0°C for 24 h. Alternatively, the cells can be mechanically disrupted using a glass/Teflon tissue grinder and allowed to extract at 0°C for 24 h. If after disrupting the cells, it is necessary to rinse the tissue grinder, or mortar and pestle, then a known volume of 90 % acetone, measured using a Class A volumetric pipette, should be used. The ease with which the pigments are removed from the cells varies considerably with different phytoplankton. In all cases, freezing the sample filters in liquid nitrogen improves extraction efficiency.

Prior to analysis, pigment extracts are vortexed and centrifuged to minimize cellular debris. To remove fine glass fiber and cellular debris from the extract, as well as enhance the life expectancy of the HPLC column, filter the extract through 13 mm PTFE (polytetrafluoroethylene) membrane syringe filters (0.2  $\mu\text{m}$  pore size). The use of Nylon filters is not recommended as they may bind certain hydrophobic pigments.

### Apparatus

The HPLC system consists of solvent pumps, sample injector, guard and analytical columns, absorption (and fluorescence) detector, and a computer. A temperature-controlled autosampler is optional, but highly recommended, to chill the samples chilled prior to injection and to reduce uncertainties during sample preparation and injection. A variety of companies manufacture HPLC systems (*e.g.* Agilent Technologies, Beckman, ThermoQuest, Waters Associates). For a review of hardware and software requirements for measuring chlorophylls and their degradation products, as well as carotenoids, see Jeffrey *et al.* (1997).

### HPLC Eluants and Gradient Programs

There are several currently recognized HPLC methods for separating chlorophylls, chlorophyll derivatives and taxonomically important carotenoids. The C<sub>18</sub> method of Wright *et al.* (1991) is recommended by SCOR and separates more than 50 chlorophylls, carotenoids, and their derivatives using a ternary gradient system. This HPLC method is described in detail below. The separation of the various pigments requires about 30 minutes. Prior to injection, 1000 µL of the aqueous acetone pigment extract is diluted with 300 µL HPLC-grade water to increase the affinity of pigments for the column during the loading step. This procedure results in sharper peaks, allowing greater loading than can be obtained with undiluted samples.

This method does not separate monovinyl and divinyl chlorophylls *a* and *b*. The presence of divinyl chlorophylls *a* and *b*, can cause errors if they are not separated either physically on the column, or by a channels ratio method from the monovinyl forms. Latasa *et al.* (1996) showed that the use of a single response factor (only for monovinyl chlorophyll *a*) could result in a 15 % to 25 % overestimation of total chlorophyll *a* concentration if divinyl chlorophyll *a* was present in significant concentrations. Although monovinyl and divinyl chlorophyll *a* co-elute, each compound absorbs differently at 436 nm and 450 nm and it is therefore possible to deconvolve the absorption signals due to these pigments (Latasa *et al.* 1996).

Alternatively, these two chlorophyll species can be separated chromatographically and individually quantified using the C<sub>8</sub> HPLC techniques described by Goericke and Repeta (1993) and Van Heukelem and Thomas (2001). The latter technique uses a two solvent system and elevated column temperature to achieve desired separations.

Regardless of the method or column-packing material used (C<sub>18</sub> or C<sub>8</sub>), it is important that HPLC performance be validated before and during use. This would include validation that resolution between peaks is acceptable, or when peaks are not chromatographically resolved, that equations based on spectral deconvolution are possible in order to quantify relative proportions of each pigment in a co-eluting pair.

### Determination of Algal Chlorophyll and Carotenoid Pigments by HPLC (Wright *et al.* 1991):

#### a. Equipment and reagents:

1. *Reagents*: HPLC grade acetone (for pigment extraction); HPLC-grade water, methanol, acetonitrile and ethyl acetate; 0.5 M ammonium acetate aq. (pH = 7.2); and BHT (2,6-di-tert-butyl-p-cresol, Sigma Chemical Co.).
2. *High-pressure injector valve* equipped with a 200µL sample loop.
3. *Guard-column* (50 mm x 4.6 mm, ODS-2 Spherisorb C<sub>18</sub> packing material, 5 µm particle size) for extending the life of the primary column.
4. *Reverse-phase HPLC column* with end capping (250 mm x 4.6 mm, 5 µm particle size, ODS-2 Spherisorb C<sub>18</sub> column).
5. *Variable wavelength or filter absorbance detector* with low volume flow through cell. Detection wavelengths are 436 nm and 450 nm.
6. *Data recording device*: a strip chart recorder, or preferably, an electronic integrator and computer equipped with hardware and software for chromatographic data analysis.
7. *Glass syringe (500 µL) or HPLC autosampler*.

8. *HPLC Solvents*: solvent A (80:20, by volume; methanol:0.5 M ammonium acetate aq., pH=7.2; 0.01 % BHT, w:v), solvent B (87.5:12.5, by volume; acetonitrile:water; 0.01 % BHT, w:v) and solvent C (ethyl acetate). Solvents A and B contain BHT to prevent the formation of chlorophyll *a* allomers. Use HPLC-grade solvents. Measure volumes before mixing. Filter solvents through a solvent resistant 0.4  $\mu\text{m}$  filter before use, and degas with helium, or an in-line vacuum degassing system, during analysis.
9. *Calibration standards*: Chlorophylls *a* and *b* and  $\beta$ , and  $\beta$ -carotene can be purchased from Sigma Chemical Co. (St. Louis, MO 63178, USA). Other pigment standards can be purchased from the International Agency for  $^{14}\text{C}$  Determination, VKI Water Quality Institute, Agern Allé 11, DK-2970 HØrsholm, Denmark. The concentrations of all standards in the appropriate solvents should be determined, using a monochromator-based spectrophotometer, prior to calibration of the HPLC system (Latasa *et al.* 1999). Spectrophotometric readings should be made at a bandwidth  $\leq 2$  nm and the optical density (*OD*) of the pigment standards should range between 0.2 to 0.8 OD units at  $\lambda_{\text{max}}$  (Marker *et al.* 1980). The recommended extinction coefficients for the various phytoplankton pigments can be found in Appendix E of Jeffrey *et al.* (1997). Absorbance is measured in a 1 cm cuvette at the peak wavelength  $\lambda_{\text{max}}$ , and at 750 nm to correct for light scattering.
10. Concentrations of the standards are calculated as

$$C_{\text{STD}}^i = \frac{10^6 [A^i(\lambda_{\text{max}}^i) - A^i(750)]}{bE_{1\text{cm}}^i}, \quad (2.1)$$

where  $C_{\text{STD}}^i$  is the concentration ( $\mu\text{g L}^{-1}$ ) of the standard for pigment *i*,  $A^i(\lambda_{\text{max}}^i)$  and  $A^i(750)$  are absorbances at  $\lambda_{\text{max}}^i$  and 750 nm, respectively, *b* is the pathlength of the cuvette (cm), and  $E_{1\text{cm}}^i$  is the weight-specific absorption coefficient ( $\text{L g}^{-1} \text{cm}^{-1}$ ) of pigment *i*. Values for  $\lambda_{\text{max}}^i$  and  $E_{1\text{cm}}^i$  are given in Appendix E of Jeffrey *et al.* (1997). Standards stored under nitrogen in the dark at  $-20^\circ\text{C}$  do not change appreciably over a one-month period, provided that they are stored in containers proven to prevent evaporation (*e.g.* glass or Teflon bottles/vials).

#### b. Procedure:

1. Set up and equilibrate the HPLC system with eluant A at a flow rate of  $1 \text{ mL min}^{-1}$ .
2. Calibrate the HPLC system using working standards prepared, on the day of use, by diluting the primary standard with the appropriate solvent (Jeffrey *et al.* 1997, Appendix E). When preparing calibration standards, one should only use dilution devices for which the precision and uncertainty have been validated with the solvent to be measured. Prepare at least 5 concentrations ( $\mu\text{g L}^{-1}$ ) of working standards for each pigment spanning the concentration range appropriate for the samples to be analyzed.
3. For each working standard, mix 1000  $\mu\text{L}$  with 300  $\mu\text{L}$  of distilled water, shake, and equilibrate for 5 min prior to injection (diluting the standards and sample extracts with water increases the affinity of pigments for the column in the loading step, resulting in an improved separation of the more polar pigments). Rinse the sample syringe twice with 300  $\mu\text{L}$  of the diluted working standard and draw 500  $\mu\text{L}$  of the working standard into the syringe for injection. Place the syringe in the injector valve, overfilling the 200  $\mu\text{L}$  sample loop 2.5-fold. To check for possible interferences in the extraction solvent and/or filter, prepare a blank by extracting a glass fiber filter in 90 % acetone, mixing 1000  $\mu\text{L}$  of the 90 % acetone filter extract and 300  $\mu\text{L}$  distilled water, and injecting the mixture onto the HPLC system. For each pigment *i*, plot absorbance peak areas (arbitrary system units) against working standard pigment masses (concentrations multiplied by injection volume). The HPLC system response factor  $F^i$  ( $\text{area } \mu\text{g}^{-1}$ ) for pigment *i* is calculated as the slope of the regression of the peak areas of the parent pigment (plus areas of peaks for structurally-related isomers if present) against the pigment masses of the injected working standards ( $\mu\text{g}$ ). Structurally related isomers (*e.g.* chlorophyll *a* allomer) contribute to the absorption signal of the standards and disregarding them will result in the over-estimation of analytes in sample extracts (Bidigare 1991).

4. Prepare pigment samples for injection by mixing a 1000  $\mu\text{L}$  portion of the aqueous acetone pigment extract and 300  $\mu\text{L}$  distilled water, shake, and equilibrate for 5 min prior to injection. Inject the sample onto the HPLC column. Samples that are pre-mixed with distilled water (or other injection buffer) should not be allowed to reside in autosampler compartments for extended durations, because hydrophobic pigments will precipitate out of solution (Mantoura *et al.* 1997). For additional information regarding HPLC method implementation and injection conditions see Wright and Mantoura (1997).
5. Following injection of the sample onto the HPLC system, use the following solvent system program to separate the chlorophyll and carotenoid pigments: 0.0' (90%A, 10%B), 1.0' (100%B), 11.0' (78%B, 22%C), 27.5' (10%B, 90%C), 29.0' (100%B), and 30.0' (100%B). Degas solvents with helium or an in-line vacuum degassing system during analysis. It should be noted that method performance varies significantly between HPLC systems because of differences in dwell volume, equilibration time, and injection conditions. It is, therefore, recommended that analysts validate that desired peak separations are attained for pigment pairs of interest by calculating the peak resolution indices  $R_s$  as

$$R_s = \frac{2(t_{R2} - t_{R1})}{w_{B1} + w_{B2}}, \quad (2.2)$$

where  $t_{R1}$  and  $t_{R2}$  are the retention times (min) of peaks 1 and 2, and  $w_{B1}$  and  $w_{B2}$  are the widths (min) of peaks 1 and 2 at their respective bases (Wright 1997). Peak separation values  $R_s < 1.0$  are insufficient for accurate quantification of peak areas (Wright 1997).

6. Peak identities are routinely determined by comparing the retention times of sample peaks with those of pure standards. Peak identities can be confirmed spectrophotometrically by collecting eluting peaks from the column outlet (or directly with an on-line diode array spectrophotometer). Absorption maxima for the various phytoplankton pigments can be found in Part IV of Jeffrey *et al.* (1997).
7. Calculate individual pigment concentrations as

$$C_{\text{Sample}}^i = \frac{A_{\text{Sample}}^i V_{\text{Extracted}} A_{\text{STD}}^{\text{Cantha}}}{F^i V_{\text{Injected}} V_{\text{Sample}} A_{\text{Sample}}^{\text{Cantha}}}, \quad (2.3)$$

where  $C_{\text{Sample}}^i$  is the individual pigment concentration ( $\mu\text{g L}^{-1}$ ),  $A_{\text{Sample}}^i$  is the area of individual pigment peak for a sample injection,  $V_{\text{Extracted}}$  is the volume extracted (mL, to nearest 0.01 mL)<sup>2</sup>,  $V_{\text{Injected}}$  is the volume injected (mL, measured to the nearest 0.001 mL),  $V_{\text{Sample}}$  is the sample volume filtered (L, measured to the nearest 0.001 L), and the other coefficients are defined above.

8. This method is designed for the separation of chlorophyll and carotenoid pigments, but it is also capable of separating the major chlorophyll breakdown products.
9. The uncertainty of the HPLC method was assessed by performing triplicate injections of a mixture of phytoplankton and plant extracts; coefficients of variation (standard deviation/mean x 100 %) ranged from 0.6 % to 6.0 %. The use of an appropriate internal standard, such as canthaxanthin, will decrease the uncertainty.

## 2.4 QUALITY ASSURANCE PROCEDURES

Quality assurance procedures outlined here should be routinely employed to insure accurate, precise and representative results.

<sup>2</sup> If an internal standard, such as canthaxanthin, is not used in the HPLC analysis, then 0.22 mL (Trees 1978; or a value determined independently by the analyst) should be added to  $V_{\text{Extracted}}$  to account for water retention by the GF/F filter (Section 2.3 above).

As a means of monitoring an instrument's performance, individual pigment response factors ( $F^i$ ) should be charted as functions of time (Clesceri *et al.* 1998). These quality control graphs should be retained with the data analysis logbooks to document the quality of each data set.

A selected number of samples should be analyzed in duplicate (or triplicate) to assess representativeness and uncertainty in the method and instrumentation. In multi-ship/investigator studies, replicate samples should be collected and archived for future intercalibration checks.

Fortified samples should be analyzed as part of the quality assurance effort. Fortified samples are prepared in duplicate by spiking a sample with known quantities of the analytes of interest at concentrations within the range expected in the samples. Fortified samples are used to assess the method's uncertainty in the presence of a typical sample matrix.

The method detection limit (*MDL*) for the analytes of interest can be determined by measuring seven replicate standard injections (Glaser *et al.* 1981). The standard deviation  $S_c$  of the seven replicate measurements is calculated, and the *MDL* is computed as

$$MDL = t(6, 0.99)S_c. \quad (2.4)$$

where  $t(6, 0.99)$  is the Student's *t* value for a one-tailed test at the 99 % confidence level, with (N-1)=6 degrees of freedom. For this particular sample size (N=7) and the 99% confidence level,  $t(6, 0.99) = 3.707$  (Abramowitz and Segun 1968, Table 26.10).

System and spiked blanks should be routinely analyzed. A system blank consists of a filter, reagents, and the glassware and hardware utilized in the analytical scheme. The system blank is quantified under identical instrumental conditions as the samples and is analyzed by appropriate quantitative methods. The system blank may not contain any of the analytes of interest above the *MDL* or corrective action must be taken. A spiked blank is defined as a system blank plus an authentic external standard containing the analytes of interest. Each set of samples should be accompanied by a spiked blank and is quantified under the same instrumental conditions as the samples.

## 2.5 PROTOCOL STATUS AND FUTURE DIRECTIONS FOR RESEARCH

Recent studies have identified the presence of novel bacterial phototrophs in coastal and oceanic waters. These include proteorhodopsin-containing *Bacteria* (Béjà *et al.* 2000, 2001) and aerobic anoxygenic phototrophic *Bacteria* (Kolber *et al.* 2000, 2001). Sequence analysis of BAC clone libraries prepared from Monterey Bay, Station ALOHA and the Southern Ocean revealed that numerous uncultivated members of the  $\gamma$ -*Proteobacteria* contain genes that code for proteorhodopsin. This membrane-bound pigment contains *trans*-retinal, absorbs at blue-green to green wavelengths, and functions as a light-driven proton pump. In an unrelated study, Kolber *et al.* (2000) used an infrared fast repetition rate (IRFRR) fluorometer to document the widespread occurrence of aerobic anoxygenic phototrophs (AAPs) in the world oceans. These microbes possess low amounts of bacteriochlorophyll *a* ( $\lambda_{max} = 358$  nm, 581 nm and 771 nm) and unusually high levels of bacteriocarotenoids ( $\lambda_{max} = 454$  nm, 465 nm, 482 nm and 514 nm). They require molecular oxygen for growth. One of us (RRB) has initiated HPLC pigment analysis of these latter clones and retinal-related compounds to determine if the Wright *et al.* (1991) method can be used for their separation and quantification.

## REFERENCES

- Abramowitz, A. and I.A. Segun, 1968: Handbook of Mathematical Functions, Dover, New York (5<sup>th</sup> Printing), 1046pp.
- Andersen, R.A., R.R. Bidigare, M.D. Keller, and M. Latasa, 1996: A comparison of HPLC pigment signatures and electron microscopic observations for oligotrophic waters of the North Atlantic and Pacific Oceans. *Deep-Sea Res. II*, **43**, 517-537.
- Béjà O, L. Aravind, E. V. Koonin, M. T. Suzuki, A. Hadd, L. P. Nguyen, S. B. Jovanovich, C. M. Gates, R. A. Feldman, J. L. Spudich, E. N. Spudich, and E. F. DeLong, 2000: Bacterial rhodopsin: Evidence for a new type of phototrophy in the sea. *Science*, **289**, 1902-1906.

- Béjà, O., E. N. Spudich, J. L. Spudich, M. LeClerc, and E. F. DeLong, 2001: Proteorhodopsin phototrophy in the ocean. *Nature*, **411**, 786-789.
- Bianchi, T. S., C. Lambert, and D. C. Biggs. 1995: Distribution of chlorophyll *a* and pheopigments in the northwestern Gulf of Mexico: a comparison between fluorometric and high-performance liquid chromatography measurements. *Bull. Mar. Science* **56**, 25-32.
- Bidigare, R.R., 1991: Analysis of algal chlorophylls and carotenoids. In: *Marine Particles: Analysis and Characterization*, D.C. Hurd and D.W. Spencer, Eds., Am. Geophys. Union, Washington, DC, 119-123.
- Bidigare, R.R., and M.E. Ondrusek, 1996: Spatial and temporal variability of phytoplankton pigment distributions in the central equatorial Pacific Ocean. *Deep-Sea Res. II*, **43**, 809-833.
- Brock, T.D., 1983: *Membrane filtration: a user's guide and reference manual*. Science Tech., Madison, WI, 381 pp.
- Chavez, F., K.R. Buck, R.R. Bidigare, D.M. Karl, D. Hebel, M. Latasa, L. Campbell, and J. Newton, 1995: On the chlorophyll *a* retention properties of glass-fiber GF/F filters. *Limnol. Oceanogr.*, **40**, 428-433.
- Clesceri, L.S., A.E. Greenberg, and A.D. Eaton (editors), 1998: Part 10000, Biological Examination, Section 1020 B. in *Standard Methods for the Examination of Water and Wastewater*. 20<sup>th</sup> ed. Baltimore (MD): American Public Health Association, American Water Works Association, Water Environment Federation.
- Dickson, M.-L., and P.A. Wheller, 1993: Chlorophyll *a* concentrations in the North Pacific: Does a latitudinal gradient exist? *Limnol. Oceanogr.*, **38**, 1813-1818.
- Gibb, S.W., R.G. Barlow, D.G. Cummings, N.W. Rees, C.C. Trees, P. Holligan and D. Suggett, 2000: Surface phytoplankton pigment distribution in the Atlantic: an assessment of basin scale variability between 50°N and 50°S. *Progress in Oceanography*, **45** (3-4), 329-368.
- Glaser, J.A., D.L. Foerst, G.D. McKee, S.A. Quave, and W.L. Budde, 1981: Trace analyses for wastewaters. *Environ. Sci. Technol.*, **15**, 1426-1435.
- Goericke, R., and D.J. Repeta, 1993: Chlorophylls *a* and *b* and divinyl chlorophylls *a* and *b* in the open subtropical North Atlantic Ocean. *Mar. Ecol. Prog. Ser.*, **101**, 307-313.
- Gordon, H.R., and D.K. Clark, 1980: Remote sensing optical properties of a stratified ocean: an improved interpretation. *Appl. Optics*, **19**, 3,428-3,430.
- Hoepffner, N., and S. Sathyendranath, 1992: Bio-optical characteristics of coastal waters: absorption spectra of phytoplankton and pigment distribution in the western North Atlantic. *Limnol. Oceanogr.* **37**, 1660-1679.
- Holm-Hansen, O., C.J. Lorenzen, R.W. Holmes, and J.D.H. Strickland, 1965: Fluorometric determination of chlorophyll. *J. du Cons. Intl. Pour l'Expl. de la Mer.*, **30**, 3-15.
- Jeffrey, S.W., and G.F. Humphrey, 1975: New spectrophotometric equations for determining chlorophylls *a*, *b*, *c*<sub>1</sub> and *c*<sub>2</sub> in higher plants, algae and natural phytoplankton. *Biochem. Physiol. Pflanzen*, **167**, 191-194.
- Jeffrey, S.W., R.F.C. Mantoura, and S.W. Wright (eds.), 1997: *Phytoplankton Pigments in Oceanography*, Monographs on Oceanographic Methodology, UNESCO, 661 pp.
- Kolber, Z. S., C. L. Van Dover, R. A. Niederman, and P. G. Falkowski, 2000: Bacterial photosynthesis in surface waters of the open ocean. *Nature*, **407**, 177-179.
- Kolber, Z. S., F. G. Plumley, A. S. Lang, J. T. Beatty, R. E. Blankenship, C. L. VanDover, C. Vetriani, M. Koblizek, C. Rathgeber, and P. G. Falkowski, 2001: Contribution of aerobic photoheterotrophic bacteria to the carbon cycle in the ocean. *Science*, **292**, 2492-2495.
- Latasa, M., R. R. Bidigare, M. E. Ondrusek, and M. C. Kennicutt II, 1996: HPLC analysis of algal pigments: A comparison exercise among laboratories and recommendations for improved analytical performance. *Mar. Chem.*, **51**, 315-324.
- Latasa, M., R. R. Bidigare, M. E. Ondrusek, and M. C. Kennicutt II, 1999: On the measurement of pigment concentrations by monochromator and diode-array spectrophotometers. *Mar. Chem.*, **66**, 253-254.

- Letelier, R.M., R.R. Bidigare, D.V. Hebel, M.E. Ondrusek, C.D. Winn, and D.M. Karl, 1993: Temporal variability of phytoplankton community structure at the U.S.-JGOFS time-series Station ALOHA (22°45'N, 158°W) based on HPLC pigment analysis. *Limnol. Oceanogr.*, **38**, 1,420-1,437.
- Mantoura, R.F.C., R.G. Barlow and E.J.H. Head, 1997: Simple isocratic HPLC methods for chlorophylls and their degradation products. Ch. 11 in Jeffrey, S.W., R.F.C. Mantoura, and S.W. Wright (editors), *Phytoplankton pigment in oceanography: guidelines to modern methods*. Vol. 10, Monographs on oceanographic methodology. UNESCO Publishing, 661 pp.
- Marker, A.F.H., E.A. Nusch, H. Rai and B. Riemann, 1980: The measurement of photosynthetic pigments in freshwaters and standardization of methods: conclusion and recommendations. *Arch. Hydrobiol. Beih. Ergebn. Limnol.* **14**: 91-106.
- Phinney, D.A., C.S. Yentsch, 1985: A novel phytoplankton chlorophyll technique: Toward automated analysis. *J. Plankton Res.*, **7**, 633-642.
- Smith, R. C., R. R. Bidigare, B. B. Prezelin, K. S. Baker, and J. M. Brooks, 1987: Optical characterization of primary productivity across a coastal front. *Mar. Biol.* **96**, 575-591.
- Snyder, L.R. and Kirkland, J.J., 1979: Quantitative and trace analysis. In: *Introduction to modern liquid chromatography*, John Wiley and Sons, New York, 541-574.
- Strickland, J.D.H., and T.R. Parsons, 1972: *A Practical Handbook of Sea Water Analysis*, Fisheries Research Board of Canada, 310 pp.
- Tester, P. A., M. E. Geesey, C. Guo, H. W. Paerl, and D. F. Millie, 1995: Evaluating phytoplankton dynamics in the Newport River estuary (North Caroline, USA) by HPLC-derived pigment profiles. *Mar. Ecol. Prog. Ser.* **124**, 237-245.
- Trees, C.C., 1978: Analytical analysis of the effect of dissolved solids on suspended solids determination. *J. Water Pollution Control Federation*, Oct 1978. pp 2370-2373.
- Trees, C.C., M.C. Kennicutt II, and J.M. Brooks, 1985: Errors associated with the standard fluorometric determination of chlorophylls and pheopigments. *Mar. Chem.*, **17**, 1-12.
- Trees, C.C., D.C. Clark, R.R. Bidigare, M.E. Ondrusek and J.L. Mueller, 2000. Accessory pigments versus chlorophyll *a* concentrations within the euphotic zone: a ubiquitous relationship. *Limnol. Oceanogr.*, **45**(5): 1130-1143.
- UNESCO, 1994: Protocols for the Joint Global Ocean Flux Study (JGOFS) Core Measurements, Manual and Guides **29**, 170pp.
- Van Heukelem, L. and C.S. Thomas, 2001: Computer-assisted high-performance liquid chromatography method development with applications to the isolation and analysis of phytoplankton pigments. *J. Chrom. A.* **910**:31-49.
- Wright, S.W., S.W. Jeffrey, R.F.C. Mantoura, C.A. Llewellyn, T. Bjornland, D. Repeta, and N. Welschmeyer, 1991: Improved HPLC method for the analysis of chlorophylls and carotenoids from marine phytoplankton. *Mar. Ecol. Prog. Ser.*, **77**, 183-196.
- Wright, S.W., 1997: Summary of terms and equations used to evaluate HPLC chromatograms. Appendix H in Jeffrey, S.W., R.F.C. Mantoura, and S.W. Wright (editors), *Phytoplankton pigment in oceanography: guidelines to modern methods*. Vol. **10**, Monographs on oceanographic methodology. UNESCO Publishing, 661 pp.
- Wright, S.W., and R.F.C. Mantoura, 1997: Guidelines for selecting and setting up an HPLC system and laboratory. Ch. 15 in Jeffrey, S.W., R.F.C. Mantoura, and S.W. Wright (editors), *Phytoplankton pigment in oceanography: guidelines to modern methods*. Vol. **10**, Monographs on oceanographic methodology. UNESCO Publishing, 661 pp.

## Chapter 3

# Fluorometric Chlorophyll *a*: Sampling, Laboratory Methods, and Data Analysis Protocols

Charles C. Trees<sup>1</sup>, Robert R. Bidigare<sup>2</sup>, David M. Karl<sup>2</sup> Laurie Van Heukelem<sup>3</sup>  
and John Dore<sup>2</sup>

<sup>1</sup>*Center for Hydro-Optics & Remote Sensing, San Diego State University, California*

<sup>2</sup>*Department of Oceanography, University of Hawaii, Hawaii*

<sup>3</sup>*Horn Point Laboratory, University of Maryland Center for Environmental Science, Horn Point, Maryland*

### 3.1 INTRODUCTION

In addition to HPLC analyses, it is recommended that the standard fluorometric methodology used for measuring chlorophylls and pheopigments also be applied to (i) the same extracted pigment samples used for HPLC analysis, and (ii) additional independent samples. Analysis of fluorometric chlorophyll *a* concentration is a far simpler procedure than HPLC analysis, especially at sea. On a given research cruise, therefore, it is economically feasible to acquire and process many more fluorometric than HPLC samples and to statistically relate fluorometric and HPLC chlorophyll *a* concentrations using linear regression analysis. This additional analysis will also enable a direct link to the historical bio-optical algorithms and database development during the CZCS validation experiments.

Protocols for fluorometric determination of the concentrations of chlorophyll and pheopigments were developed initially by Yentsch and Menzel (1963) and Holm-Hansen *et al.* (1965), and are described in detail by Strickland and Parsons (1972). Holm-Hansen *et al.* (1965) and Strickland and Parsons (1972) used first principles of fluorescence spectroscopy to derive these fluorometric equations. The equation proposed by Yentsch and Menzel (1963) is only indirectly linked to first principles, through debatable assumptions, and its use is not recommended. Although these measurements have been shown to contain errors as compared to HPLC determinations (Trees *et al.* 1985; Smith *et al.* 1987; Hoepffner and Sathyendranath 1992; Bianchi *et al.* 1995; Tester *et al.* 1995), the CZCS phytoplankton pigment concentration algorithms were based on them entirely. The SeaWiFS protocols for this analysis will be those given in Strickland and Parsons (1972) as updated by this chapter.

Pigment databases generally show a log-normal distribution, which is consistent with that proposed by Campbell (1995) for bio-optical properties. Therefore, it is appropriate to perform log-linear regressions on HPLC determined total chlorophyll *a* (chlorophyllide *a*, chlorophyll *a* epimer, chlorophyll *a* allomer, monovinyl chlorophyll *a* and divinyl chlorophyll *a*) and fluorometrically determined chlorophyll *a*, using model I regressions. Standard Model I regressions were selected because HPLC determined total chlorophyll *a* concentrations are to be predicted from fluorometrically determined chlorophyll [Model I regressions are appropriate for both predictions and determining functional relationships, whereas Model II regressions should not be used to predict values of *y* given *x* (page 543, Sokal and Rohlf 1995)].

Examples of regression models predicting log HPLC total chlorophyll *a* (following Chapter 2 HPLC protocols) from log fluorometric chlorophyll *a* are shown in Figures 3.1, 3.2, and 3.3 for three cruises in different geographic areas. In each example, the regression slopes are significantly different from a one-to-one relationship, although for the Gulf of California (GoCAL November 1996, Figure 3.3) the slope is close to unity. One-to-one ratios have also been found for other geographic areas, but not necessarily during all seasons. Therefore, the relationship (slope and offset) between HPLC total chlorophyll *a* and fluorometric chlorophyll *a* must be determined for a selected number of samples for each cruise, so that a cruise-specific scaling factor can be applied to other fluorometric samples.

The protocols specified below for fluorometric chlorophyll *a* analyses follow closely those prescribed in the *JGOFS Core Measurement Protocols* (UNESCO 1994), but they differ in one important respect. Absorption of light in seawater, or any other medium, is a volumetric process, even though the volume absorption coefficient may

vary with the density of the medium. For ocean color and optical analyses, therefore, the concentration of chlorophyll *a* shall be expressed in units of mass per unit volume of seawater, either in  $\mu\text{g L}^{-1}$ , or  $\text{mg m}^{-3}$ . This differs from the JGOFS protocols, which specify that concentrations in seawater of chlorophyll *a* and pheopigments should be expressed in  $\mu\text{g kg}^{-1}$ .

### 3.2 SAMPLE ACQUISITION AND STORAGE

Water samples should be taken using, *e.g.*, Niskin bottles at the site of, and simultaneously with, the surface in-water upwelled radiance and reflectance measurements, and at depth increments sufficient to resolve variability within at least the top optical depth.

The  $K(z)$ , profiles over this layer will be used to compute optically weighted, near-surface pigment concentration for bio-optical algorithm development (Gordon and Clark 1980). When possible, samples should also be acquired at several depths distributed throughout the upper 200 m of the water column [or in turbid water, up to seven diffuse attenuation depths, *i.e.*  $\ln(E(0)/E(z))=7$ , to provide a basis for relating fluorescence signals to pigment mass concentration.

Samples should be filtered as soon as possible after collection. If processing must be delayed for more than an hour, hold the samples on ice, or in a freezer at 4°C, and protect them from exposure to light. For delays longer than several hours, the samples should be stored in liquid nitrogen. Use opaque sample bottles, because even brief exposure to light during sampling and/or storage might alter pigment values.

#### *Filtration*

Whatman GF/F glass fiber filters, with approximately 0.7  $\mu\text{m}$  pore size, are preferred for removing phytoplankton from water. The glass fibers assist in breaking the cells during grinding and no precipitate forms after acidification. Twenty-five mm diameter GF/F glass fiber filters should be used with a vacuum or positive pressure with a pressure differential equivalent to 180-200 mm of mercury. Large filtration volumes are not required, because of the increased sensitivity of the fluorescence measurement.

Inert membrane filters, such as polyester filters, may be used when size fraction filtration is required. When this is done, it is recommended to also filter a replicate sample through a GF/F to determine the total concentration. Summing the various size-fractionated concentrations may not produce an accurate estimate of the total, because of the potential for cell disruption during filtration.

There has been an ongoing discussion on filter types and retention efficiencies for natural samples. Phinney & Yentsch (1985) showed the inadequacy of GF/F filters for retaining chlorophyll *a* in oligotrophic waters, as did Dickson and Wheeler (1993) for samples from the North Pacific. In response to Dickson and Wheeler (1993), Chavez *et al.* (1995) compared samples collected in the Pacific Ocean using GF/F and 0.2  $\mu\text{m}$  membrane filters with small filtered volumes (100-540 mL). Their results for small volumes showed a very close agreement between the two filter types with GF/F filters having only a slightly positive 5% bias.

Filtration volume can directly affect the retention efficiency for GF/F filters. Particles can be retained by filters through a variety of ways, such as filter sieving, filter adsorption, electrostatic and van der Waals attractions (Brock, 1983). When water flows through the pores of a Nuclepore filter, streamlines are formed that can align small particles longitudinally, with the result that cell diameter becomes important with these filters. It is known, on the other hand, that Whatman GF/F filters can retain particles much smaller than their rated pore size. Generally, at small volumes (100-300 mL) filter adsorption, and electrostatic and van der Waals attractions are important, whereas at larger volumes (> 2,000 mL) sieving dominates. This has been tested in oligotrophic waters off Hawaii in which small (< 500 mL) and large volumes (> 2-4 liters) retained similar amounts of chlorophyll *a* on the two types of filters, whereas for intermediate sample volumes the GF/F filters showed lower concentrations. As a general rule, it is recommended that the following volumes be filtered for these water types: 0.5-1.0 liter for oligotrophic, 0.2-0.5 liter for mesotrophic, and 0.1 liter and less for eutrophic water.

It is recommended to not pre-filter seawater samples to remove large zooplankton and particles, because this practice may exclude pigment-containing colonial and chain-forming phytoplankton, such as diatoms and *Trichodesmium* sp. Forceps should be used to remove large zooplankton from the GF/Fs following filtration.

### *Sample Handling, and Storage*

Samples should be filtered as quickly as possible after collection, and the filters stored immediately in liquid nitrogen. Liquid nitrogen is the best method for storing filter samples with minimum degradation for short, as well as, longer storage times (*e.g.* 1 year). Placing samples in liquid nitrogen also assists in pigment extraction by weakening the cell wall and membrane during this rapid temperature change. Ultra-cold freezers (-90°C) can be used for storage, although they have not been tested for longer than 60 days (Jeffrey *et al.* 1997). Conventional deep freezers should not be used for storing samples more than 20 hours before transferring them to an ultra-cold freezer, or liquid nitrogen.

Again, storage of samples in liquid nitrogen immediately after filtration is the preferred method. The addition of MgCO<sub>3</sub> at the end of the filtration process to stabilize chlorophyll has not been used for many years as a routine oceanographic method, because of the uncertainty in pigment absorption by MgCO<sub>3</sub>.

If samples are to be stored for any length of time prior to fluorometric analysis, they should be folded in half with the filtered halves facing in. This eliminates problems of rubbing particles off the filter during placement in sample containers and storage.

It is strongly recommended to use aluminum foil wrappings for sample containers. This simple, but effective, container is both inexpensive and easy to use. Cut small pieces of heavy-duty aluminum foil into approximately 4 cm squares. Fold each piece in half, and using a fine-point permanent marker, write a short sample identifier (*e.g.* first letter of the cruise and a sequential sample number) on the foil. Writing on the folded foil, prior to placement of the filter, both avoids puncturing the foil with the marking pen, and improves the legibility of the sample identifier. Place the folded filter in the aluminum foil. Fold the three open sides to form an envelope that is only slightly larger than the folded filter (~3cm x 1.5cm).

The use of foil containers minimizes the size requirement of the storage container. It is also acceptable to use either cryogenic tubes, or HistoPrep tissue capsules, but they occupy more storage volume per sample, and they are more expensive than aluminum foil. If fluorometric analysis is to be done soon after collection, it is still recommended to place the samples in liquid nitrogen to assist in pigment extraction, and on removal from the liquid nitrogen to place them immediately in chilled 90% acetone.

### *Recordkeeping*

Information regarding sample identification should be logged in a laboratory notebook with the analyst's initials. For each filter sample record the sample identifier (as written on the sample container), station number for the cruise, water volume filtered ( $V_{\text{FILT}}$ ) in mL, and depth of the water sample, together with the date, time, latitude, and longitude of the bottle cast during which the sample was acquired.

## **3.3 LABORATORY METHODS FOR FLUOROMETRIC DETERMINATION OF CHL. *a* AND PHEOPIGMENT CONCENTRATIONS**

Chlorophyll and pheopigments can be determined using either a Turner Designs (or Sequoia) fluorometers equipped with the standard light sources and Corning excitation and emission filters, following the manufacturer's recommendation for measuring extracted chlorophyll. The fluorometric instrument should be warmed-up for at least 30 to 45 minutes prior to making measurements.

Because of the acidification requirement for the standard fluorometric method (Holm-Hansen *et al.* 1965), differences in excitation and emission wavelength bands between fluorometers can produce uncertainties (Trees *et al.* 1985). The sensitivity with which a particular instrument is able to differentiate between chlorophyll and pheopigment is a function of the excitation wavelength. This effect is measured during calibration of the fluorometer and is called the tau factor ( $\tau$ ). Saijo and Nishizawa (1969) have shown that  $\tau$  can vary from 1 to 11.5, depending upon the excitation wavelength (in the range between 410 nm and 440 nm). For example, a comparison between a Turner Designs (Model 10-005R) analog fluorometer and a Turner Designs (Model 10-AU-005) digital fluorometer showed statistically significant differences for 42 oceanic samples (slope = 1.06), even though both were calibrated with exactly the same standards (Figure 3.4). The departure from a unit slope is attributable to differences in the excitation bands for the two fluorometers.

### Fluorometer Calibrations

Bench fluorometers used to measure concentrations of extracted chlorophyll and pheopigments should be calibrated using authentic chlorophyll *a* standards, as prescribed also in the HPLC Protocols (Chapter 2). Chlorophyll *a* standards can be purchased from Sigma Chemical Co. (St. Louis, MO 63178, USA).

If a fluorometer has been shipped for a cruise, or if it has been unused for several weeks, it is strongly recommended that it be recalibrated with an authentic chlorophyll *a* standard. The use of solid standards, like those provided by Turner Designs and other manufacturers, can only provide a check for instrumental drift. They cannot be used as primary pigment standards. However, the solid standard should be used at frequent intervals during each day's analyses to monitor instrument drift.

The concentration of the chlorophyll *a* standard, in the appropriate solvent, must be determined using a monochromator-based spectrophotometer prior to calibrating the fluorometer. The recommended extinction coefficients for chlorophyll *a* in several solvents can be found in Appendix E of Jeffrey *et al.* (1997). Absorbance is measured in a 1 cm cuvette at the peak wavelength  $\lambda_{\max}$ , and at 750 nm to correct for light scattering. The bandwidth of the spectrophotometer should be between 0.5 and 2  $\mu\text{m}$ , with the standard concentration being such that the absorbance falls between 0.1 and 1.0 optical density units (Clesceri *et al.*, 1998a). The concentration of the standard is calculated as

$$C_{\text{STD}} = \frac{10^6 [A(\lambda_{\max}) - A(750)]}{bE_{1\text{cm}}}, \quad (3.1)$$

where  $C_{\text{STD}}$  is the concentration ( $\mu\text{g L}^{-1}$ ) of the chlorophyll *a* standard,  $A(\lambda_{\max})$  and  $A(750)$  are absorbances at  $\lambda_{\max}$  and 750 nm,  $b$  is the pathlength of cuvette (cm), and  $E_{1\text{cm}}$  is the specific absorption coefficient ( $\text{L g}^{-1} \text{cm}^{-1}$ ) of chlorophyll *a* in 90% acetone. For 90% acetone  $E_{1\text{cm}} = 87.67 \text{ L g}^{-1} \text{cm}^{-1}$ , and for 100% acetone  $E_{1\text{cm}} = 88.15 \text{ L g}^{-1} \text{cm}^{-1}$ , when applied to the absorption measured at the peak wavelength  $\lambda_{\max}$  (Jeffrey *et al.* 1997, Appendix E). The peak wavelength  $\lambda_{\max}$  must be determined by inspection of the measured spectrum, because its location may shift due to interactions between the particular solvent and mixture of pigment compounds in each sample. Standards stored under nitrogen in the dark at  $-20^\circ\text{C}$  do not change appreciably over a one-month period, provided that they are stored in containers proven to prevent evaporation (*e.g.* glass or Teflon bottles/vials).

The stock chlorophyll *a* standard, with its concentration measured on a spectrophotometer as described above, should be diluted using calibrated gas-tight syringes, and Class A volumetric pipettes and flasks. The minimum number of dilutions of the stock standard for calibrating a fluorometer depends on whether it is a digital model (Turner Designs 10-AU-005), or it is an analog model with a mechanical mode for changing sensitivity (*e.g.* Turner Designs 10-005). A minimum of 5 dilutions is required for calibrating a digital fluorometer. Analog fluorometers with a variety of door settings, such as the Turner Designs Model 10-005, must be calibrated for each door setting using at least three standard concentrations per door. The diluted standard pigment concentrations used in calibrating the fluorometer must bracket the range of concentrations found in the samples being analyzed.

Each diluted chlorophyll *a* standard is placed in the fluorometer and the signal ( $F_b$ ) is recorded, after waiting a short period of time (60 seconds) for it to stabilize. The standard is removed and diluted HCL acid (2 drops of 5 %, or 1 drop of 10 %, both concentrations by volume) is added and mixed within the test tube. The tube is then placed back into the fluorometer, and after stabilization, the acidified fluorescence signal ( $F_a$ ) is recorded. Following acidification of the chlorophyll *a* standard, the fluorescence signal stabilizes relatively quickly. This is not the case for natural samples that contain a mixture of pigment compounds, however, and stabilization time may vary from sample to sample. Stabilization time has to be the same for both pigment standards and for natural samples. To minimize this source of uncertainty, and to standardize this measurement technique, it is recommended that both acidified natural sample and acidified pigment standards be allowed to react with the acid for one minute prior to recording the acidified fluorescence signal ( $F_a$ ). Two drops of 5 % (by volume) hydrochloric acid is added to each of the pigment standards and natural samples. Once the acid is added, the sample in the test tube should be mixed by inverting the tube several times, using parafilm as a stopper. All fluorometric measurements for both pigment standards and natural samples should be carried out at room temperature. A 90 % (by volume) acetone blank ( $Blk_b$ ) and an acidified acetone blank ( $Blk_a$ ) should also be measured, even though the acidified blank ( $Blk_a$ ) is frequently found to be equal to the non-acidified blank ( $Blk_b$ ). The fluorometer's sensitivity to pheopigments,  $\tau$ , is calculated as

$$\tau = \frac{F_b - Blk_b}{F_a - Blk_a}, \quad (3.2)$$

and is averaged over all concentrations of the chlorophyll *a* standard. For the mechanical door model fluorometers, data from the higher gain door settings will often become noisy and computed  $\tau$  values will begin to decrease. These data should be excluded from the average. The fluorometer's response factor,  $F_R$  ( $\mu\text{g L}^{-1}$  per fluorescence signal), is determined as the slope of the simple linear regression equation

$$C_{STD} = F_R (F_b - Blk_b), \quad (3.3)$$

calculated for the sample of diluted concentrations of the pigment standard, and forcing a zero intercept. With a digital fluorometer, the regression analysis is applied to the data from the entire 5, or more, concentrations and a single  $F_R$  factor is determined for the instrument. With a mechanical fluorometer, the regression is applied to the data from the 3, or more, concentrations of the standard, and a separate  $F_R$  factor is determined, for each door setting. As a means of monitoring an instrument's performance,  $F_R$  factors from successive calibrations should be charted as functions of time (Clesceri *et al.*, 1998b). These quality control graphs should be retained with the data analysis logbooks to document the quality of each data set for which that fluorometer is used.

#### *Solvent Preparation.*

It is recommended that 90 % acetone (by volume) be used to extract pigments for the fluorometric analysis. Richard and Thompson (1952) were the first to propose 90 % acetone as a solvent to extract pigments from marine phytoplankton. Their results indicated improved extraction efficiencies, and also showed that the procedure minimized the activity of the naturally occurring chlorophyllase enzyme, which degrades the pigment. With a graduated cylinder, make up 90 % acetone by first pouring in distilled water, followed by 100 % acetone. Using volumetric pipettes, or auto-pipettes, accurately measure 8 mL to 10 mL of 90 % acetone and place it in a centrifuge tube. Record this volume as  $V_{EXT}$ . A number of such tubes containing acetone are then stored in a freezer and individually removed as filter samples are collected. Pre-chilling the solvent in this way reduces the possibility of temperature induced pigment degradation.

#### *Extraction*

Filters are removed from liquid nitrogen and placed in the chilled centrifuge tubes for extraction in  $V_{EXT}$  mL of 90% acetone. Samples are disrupted by sonication, placed in a freezer, and allowed to extract at 0°C for 24 h. Alternatively, the cells can be mechanically disrupted using a glass/Teflon tissue grinder and allowed to extract at 0°C for 24 h. If after disrupting the cells, it is necessary to rinse the tissue grinder, or mortar and pestle, then a known volume of 90% acetone, measured using a Class A volumetric pipette, should be used. The ease at which the pigments are removed from the cells varies considerably with different phytoplankton. In all cases, freezing the sample filters in liquid nitrogen improves extraction efficiency. Prior to analysis, pigment extracts are swirled into a vortex to remove particles from the sides of the tube, and then centrifuged to minimize cellular debris.

#### *Measurement*

Following the same measurement procedure described above under *Fluorometer Calibration*, each extracted sample is placed in the fluorometer and its non-acidified and acidified responses,  $F_b$  and  $F_a$ , are measured and recorded. The concentration of chlorophyll [Chl] ( $\mu\text{g L}^{-1}$ ) in the sample is calculated as

$$[Chl] = (F_b - F_a - Blk_b + Blk_a) \frac{\tau}{\tau - 1} F_R \frac{V_{EXT}}{V_{FILT}}, \quad (3.4)$$

and pheopigments concentration [Pheo] ( $\mu\text{g L}^{-1}$ ) as

$$[Pheo] = \left\{ (F_a - Blk_a) \tau - (F_b - Blk_b) \right\} \frac{\tau}{\tau - 1} F_R \frac{V_{EXT}}{V_{FILT}}, \quad (3.5)$$

where volumes extracted  $V_{EXT}$  and filtered  $V_{FILT}$  are in mL. Pheopigment concentrations determined using the standard fluorometric method of Holm-Hansen *et al.* (1965) have not been reported in published articles for many years. This is based on the fact that (i) there is always a residual amount of pheopigments in all natural samples (Smith and Baker, 1978; 25% of the summed chlorophyll plus pheopigment), (ii) pheopigment concentrations are

overestimated in the presence of chlorophyll *b* (Lorenzen and Jeffrey, 1980; Vernet and Lorenzen, 1987), and (iii) HPLC measured pheopigments, generally contribute very little to the chlorophyll *a* pigment pool (e.g., Hallegraeff, 1981; Everitt *et al.*, 1990; and Bricaud *et al.*, 1995). Trees *et al.* (2000a) assembled an extensive HPLC pigment database (5,617 samples) extending over a decade of sampling and analysis, and including a variety of environments ranging from freshwater to marine, oligotrophic to eutrophic, and tropical to polar, and found that the average pheopigment to chlorophyll *a* ratio was only 0.037. This global scale result emphasizes the problems associated with estimating pheopigments using the standard fluorometric method.

### 3.4 *In Situ* CHLOROPHYLL *a* FLUORESCENCE PROFILES

An *in situ* fluorometer should be employed to measure a continuous profile of chlorophyll fluorescence. The fluorometer should be mounted on the same underwater package as the water sampler, ideally together with a CTD, transmissometer and other inherent optical properties (IOP) sensors. In some cases it may be desirable to also include a radiometer on this package, if shading effects associated with the package and/or ship are not significant.

*In situ* fluorometers produce nearly continuous profiles of artificially stimulated fluorescence. Fluorometer data (in volts) should be corrected by subtracting an offset, determined by shading the instrument on deck. These unscaled fluorescence responses are adequate to provide guidance in K-profile analysis and interpretation.

To produce vertical continuous profiles of pigment concentration, HPLC-derived pigment concentrations from water samples taken at discrete depths may be interpolated, with the aid of *in situ* fluorescence profiles. These *fluorescence interpolated* profiles should then be used with  $K_d(z, \lambda)$  profiles to compute the optically weighted average pigment concentration over the top attenuation length (Gordon and Clark 1980).

The A/D channel used to acquire and record signal voltages from the *in situ* fluorometer must be calibrated, and its temperature-dependent response to known voltage inputs characterized. The range dependent A/D bias coefficients should be determined at approximately 5<sup>o</sup> C intervals over the range from 0-25<sup>o</sup> C to characterize the temperature sensitivity of the data acquisition system.

Zero fluorescence offsets should be measured on deck before and after each cast; the optical windows should be shaded to avoid contamination of the zero offset value by ambient light. Before each cast, the fluorometer windows should be cleaned following the manufacturer's instructions.

### 3.5 PROTOCOL STATUS AND FUTURE DIRECTIONS FOR RESEARCH

In order to minimize interferences caused by the overlapping excitation and emission wavebands of chlorophylls *a*, *b*, *c* and pheopigments, Turner Designs (Sunnyvale, CA) manufactures the multi-spectral fluorometer TD-700. This instrument was recently tested using samples collected at the US JGOFS Hawaii Ocean Time-series Station ALOHA (22.75<sup>o</sup>N, 158<sup>o</sup>W). A set of replicate monthly (May - Dec 2000) pigment samples collected between the surface and 175 m were analyzed by HPLC using the protocols described in Chapter 2. Duplicate samples were subsequently analyzed in 100% acetone with the TD-700 using the manufacturer's calibration. The results of these comparisons are illustrated in Figures 3.5, 3.6 and 3.7 for chlorophylls *a*, *b*, and *c*, respectively. The Model I regression equations predicting each HPLC pigment (in mg m<sup>-3</sup>) from the equivalent TD700 estimate are:

- HPLC Chl *a* = 0.729[TD-700 Chl *a*] + 0.0144; ( $r^2 = 0.894$ ).
- HPLC Chl *b* = 0.607[TD-700 Chl *b*] - 0.0163; ( $r^2 = 0.816$ ).
- HPLC Chl *c* = 1.083[TD-700 Chl *c*] - 0.00249; ( $r^2 = 0.906$ ).

These equations differ significantly from a one-to-one relationship. The present comparisons differ also from those published in Trees *et al.* (2000a), although care must be used in this comparison since the concentrations were expressed there in ng L<sup>-1</sup> (which accounts for the factor of 10<sup>-3</sup> differences in the respective offset coefficients). These results call into question the stability of the fluorometer. It is also evident that the equations provided by the manufacturer must be verified with HPLC data, and that these calibration relationships should be reviewed frequently.

It is interesting and noteworthy that the TD-700 fluorometer did not detect pheopigments in any of the samples analyzed.

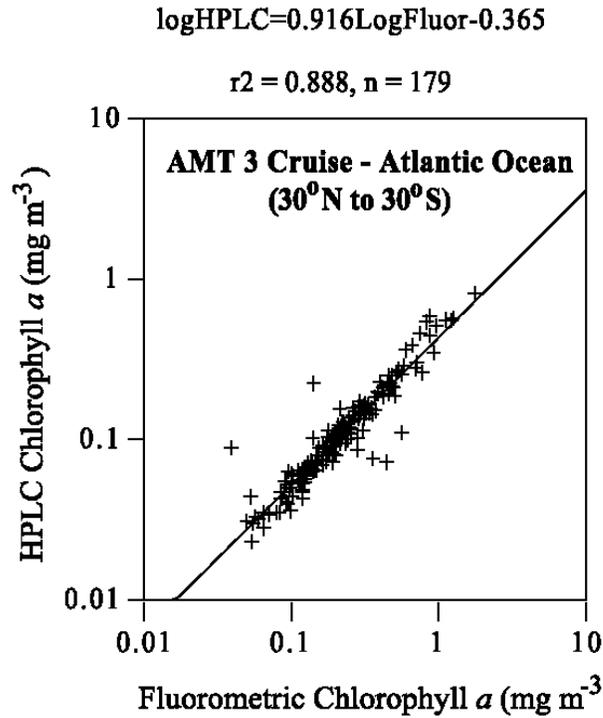


Figure 3.1: Comparisons between fluorometrically determined chlorophyll and HPLC determined total chlorophyll *a* (chlorophyllide *a*, chlorophyll *a* epimer, chlorophyll *a* allomer, monovinyl chlorophyll *a*, and divinyl chlorophyll *a*) from samples collected during Atlantic Meridional Transect 3 cruise (30°N to 30°S, October 1996).

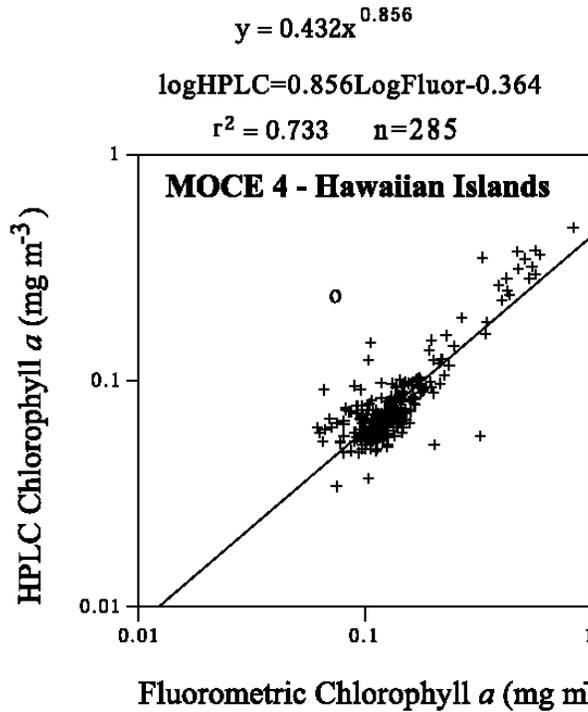


Figure 3.2: Same as Figure 3.1 for data collected during the Marine Optical Characterization Experiment (MOCE) 4 cruise.

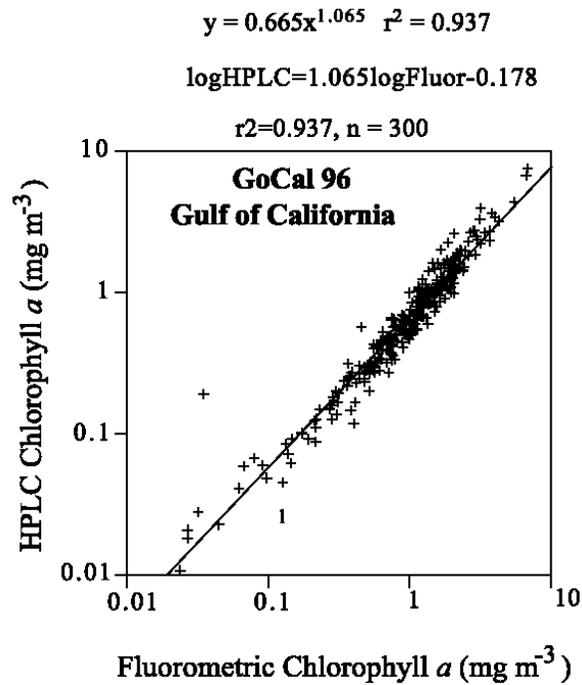


Figure 3.3 Same as Figure 3.1 for data collected during the Gulf of California cruise (Gulf of California, November 1996).

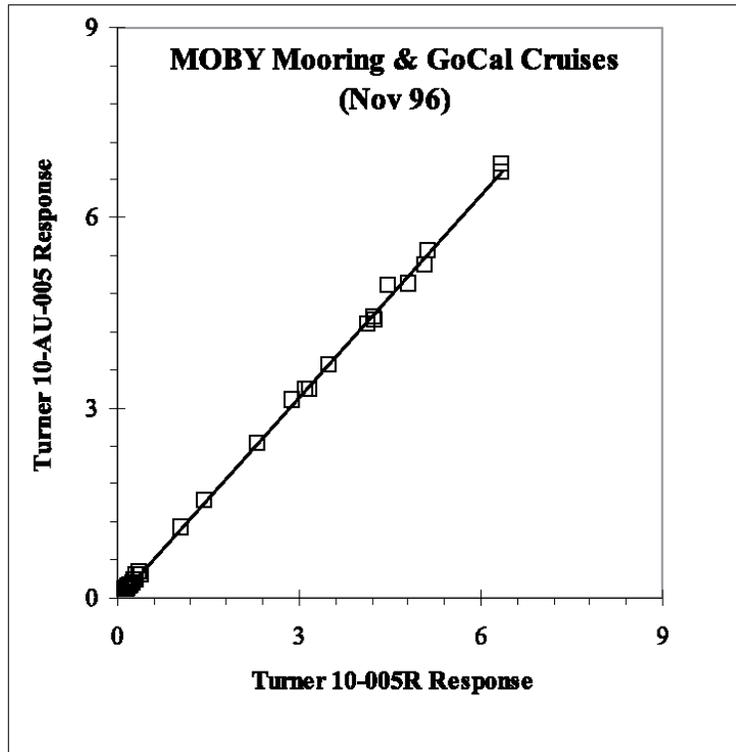


Figure 3.4. Comparison of fluorometrically determined chlorophyll *a* using the VisLab Turner Fluorometer (10-005R) and the Moss Landing Marine Laboratory Turner Fluorometer (10-AU-005). Samples were analyzed from a MOBY Nov 96 cruise and a Gulf of California cruise (Mueller, Nov 96).

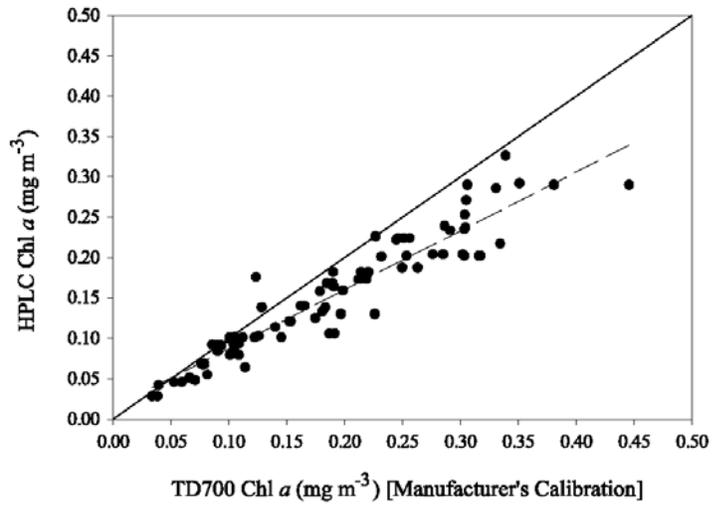


Figure 3.5. Comparison between chlorophyll a determined by the TD700 equation supplied by the manufacturer and that measured by HPLC methods.

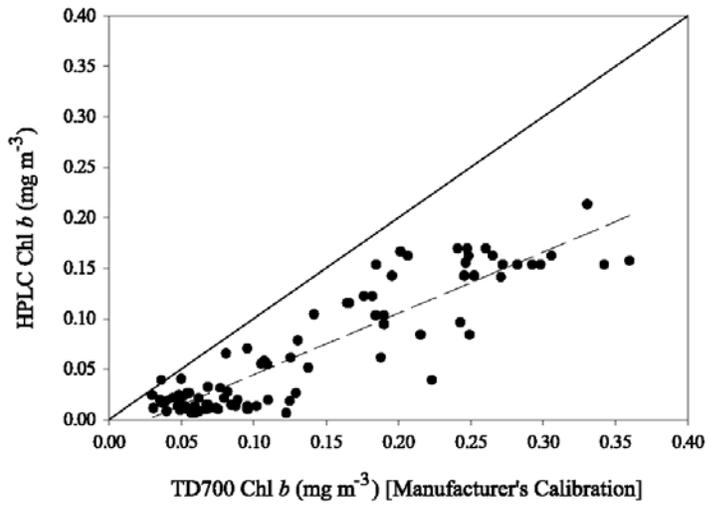


Figure 3.6. Same as Figure 3.5 for chlorophyll b.

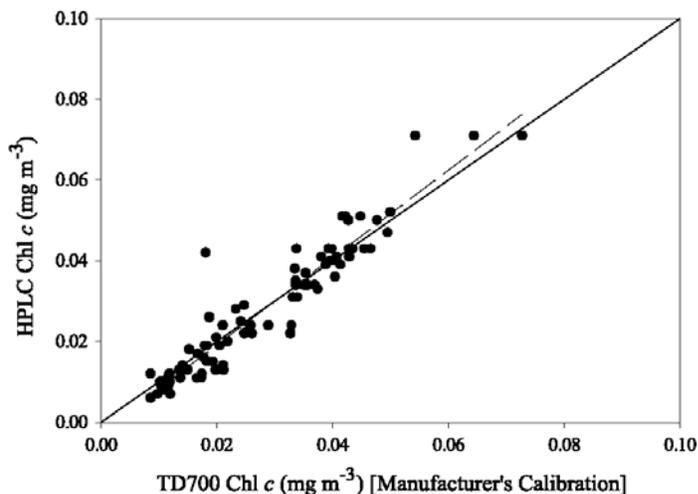


Figure 3.7. Same as Figure 3.5 for chlorophyll c.

## REFERENCES

- Bianchi, T. S., C. Lambert, and D. C. Biggs. 1995: Distribution of chlorophyll *a* and pheopigments in the northwestern Gulf of Mexico: a comparison between fluorometric and high-performance liquid chromatography measurements. *Bull. Mar. Science*, **56**, 25-32.
- Brock, T.D., 1983: *Membrane filtration: a user's guide and reference manual*. Science Tech., Madison, WI, 381 pp.
- Campbell, J.W. 1995: The lognormal distribution as a model for bio-optical variability in the sea. *J. Geophys. Res.*, **100**, 13237-13254.
- Chavez, F., K.R. Buck, R.R. Bidigare, D.M. Karl, D. Hebel, M. Latasa, L. Campbell, and J. Newton, 1995: On the chlorophyll *a* retention properties of glass-fiber GF/F filters. *Limnol. Oceanogr.*, **40**, 428-433.
- Clesceri, L.S., A.E. Greenberg and A.D. Eaton (eds), 1998a: Part 10000, Biological Examination, Section 10200 H. in *Standard Methods for the Examination of Water and Wastewater*. 20th ed. Baltimore (MD): American Public Health Association, American Water Works Association, Water Environment Federation.
- Clesceri, L.S., A.E. Greenberg and A.D. Eaton (eds), 1998b: Part 10000, Biological Examination, Section 10200 B. in *Standard Methods for the Examination of Water and Wastewater*. 20th ed. Baltimore (MD): American Public Health Association, American Water Works Association, Water Environment Federation.
- Dickson, M.-L., and P.A. Weeller, 1993: Chlorophyll *a* concentrations in the North Pacific: Does a latitudinal gradient exist? *Limnol. Oceanogr.*, **38**, 1813-1818.
- Gordon, H.R., and D.K. Clark, 1980: Remote sensing optical properties of a stratified ocean: an improved interpretation. *Appl. Optics*, **19**, 3,428--3,430.
- Hoepffner, N., and S. Sathyendranath. 1992: Bio-optical characteristics of coastal waters: absorption spectra of phytoplankton and pigment distribution in the western North Atlantic. *Limnol. Oceanogr.* **37**: 1660-1679.
- Holm-Hansen, O., C.J. Lorenzen, R.W. Holmes, and J.D.H. Strickland, 1965: Fluorometric determination of chlorophyll. *J. du Cons. Intl. Pour l'Expl. de la Mer.*, **30**, 3-15.

- Jeffrey, S.W., R.F.C. Mantoura, and S.W. Wright (eds.), 1997: *Phytoplankton Pigments in Oceanography*, Monographs on Oceanographic Methodology, UNESCO, 661 pp.
- Lorenzen, C.J. and S.W. Jeffrey. 1980: *Determination of Chlorophyll in Seawater*. UNESCO Technical Papers in Marine Science, Vol. **35**, UNESCO, 20 pp.
- Phinney, D.A. and C.S. Yentsch, 1985: A novel phytoplankton chlorophyll technique: Toward automated analysis. *J. Plankton Res.*, **7**, 633-642.
- Richards, F.A. and T.G. Thompson. 1952: The estimation and characterization of plankton populations by pigment analysis. II. A spectrophotometric method for the estimation of plankton pigments. *J. Mar. Res.*, **11**, 156-172.
- Saijo, Y. and S. Nishizawa. 1969: Excitation spectra in the fluorometric determination of chlorophyll a and phaeophytin a. *Mar Biol.*, **2**, 135-136.
- Smith, R. C., and K. S. Baker. 1978: The bio-optical state of ocean waters and remote sensing. *Limnol. Oceanogr.*, **23**, 247-259.
- Smith, R. C., R. R. Bidigare, B. B. Prezelin, K. S. Baker, and J. M. Brooks. 1987: Optical characterization of primary productivity across a coastal front. *Mar. Biol.* **96**: 575-591.
- Strickland, J.D.H., and T.R. Parsons, 1972: *A Practical Handbook of Sea Water Analysis*, Fisheries Research Board of Canada, 310 pp.
- Tester, P. A., M. E. Geesey, C. Guo, H. W. Paerl, and D. F. Millie, 1995: Evaluating phytoplankton dynamics in the Newport River estuary (North Caroline, USA) by HPLC-derived pigment profiles. *Mar. Ecol. Prog. Ser.* **124**, 237-245.
- Trees, C.C., R.R. Bidigare, D.M. Karl and L. Van Heukelem, 2000a: Fluorometric chlorophyll a: sampling, laboratory methods, and data analysis protocols, Chapter 14 in: Fargion, G.S. and J.L. Mueller (Eds.) *Ocean Optics Protocols for Satellite Ocean Color Sensor Validation*. NASA/TM-2000-209966, NASA Goddard Space Flight Center, Greenbelt, MD. pp 162-169.
- Trees, C.C., D.K. Clark, R.R. Bidigare, M.E. Ondrusek, and J.L. Mueller. 2000: Accessory pigments versus chlorophyll a concentrations within the euphotic zone: a ubiquitous relationship. *Limnol. Oceanogr.*, **45**(5): 1130-1143.
- Trees, C.C., M.C. Kennicutt II, and J.M. Brooks, 1985: Errors associated with the standard fluorometric determination of chlorophylls and pheopigments. *Mar. Chem.*, **17**, 1-12.
- UNESCO, 1994: Protocols for the Joint Global Ocean Flux Study (JGOFS) Core Measurements, Manual and Guides 29, 170pp.
- Vernet, M., and C. J. Lorenzen. 1987: The presence of chlorophyll b and the estimation of pheopigments in marine phytoplankton. *J. Plankton Res.*, **9**, 255-265.
- Yentsch, C.S., and D.W. Menzel, 1963: A method for the determination of phytoplankton, chlorophyll, and phaeophytin by fluorescence. *Deep-Sea Res.*, **10**, 221-231.



## Chapter 4

# Backscattering by Coccolithophorids and Coccoliths: Sample Preparation, Measurement and Analysis Protocols

William M. Balch and David T. Drapeau

*Bigelow Laboratory for Ocean Sciences, E. Boothbay Harbor, Maine*

### 4.1 INTRODUCTION

Estimates of backscattering are of direct relevance to understanding reflectance properties of marine phytoplankton. This is particularly true for certain phytoplankton species that are highly optically-scattering, such as coccolithophores that produce calcium carbonate scales (“coccoliths”) and shed them into seawater. Due to the high rates of coccolith production by coccolithophorids, the small size of the coccoliths (2  $\mu\text{m}$  to 5  $\mu\text{m}$ ) and low sinking rate (0.1  $\text{m d}^{-1}$ ), suspended coccoliths are ubiquitous in the world ocean, and their residence time can be considerable. They typically are responsible for 10-20% of the total backscattering in the sea, and they thus contribute a significant part of the ocean’s radiative budget. Their presence also can cause errors in the remote sensing of chlorophyll (Balch *et al.* 1989). Coccolithophores produce mesoscale blooms spanning hundreds of thousands of square kilometers, and the large bloom size makes remote sensing one of the few ways to study their spatial extent. Therefore, knowledge of coccolithophore and coccolith backscattering properties is critical for understanding their overall contributions to remote sensing reflectance, as well as for development of algorithms for detecting their abundance.

Commercially available instruments<sup>3</sup> for measuring backscattering rely on measurement of volume scattering at a single angle (141° for the HOBILABS Hydrosat; 117° for the WET Labs ECO meter) or three angles (100°, 125°, and 150°; WET Labs EcoVSF3). While one to three angles can do a reasonable job of characterizing the backward part of the volume scattering function (VSF), measurements of the VSF at more angles improves the definition of its shape, and subsequent integration over the backward directions provides a more accurate estimate of the backscattering coefficient. One limitation of *in situ* instruments is that they do not allow easy sample manipulation; a benchtop instrument allows measurements under much more controlled conditions than can be found in the field and it also allows the manipulation of samples prior to sampling (such as the experimental removal of coccoliths from the suspension). Because of the importance of calcium carbonate coccoliths to light scattering in the sea, two other ancillary measurements are important for algorithm development: microscopic counts of coccolith concentration and measurements of particulate inorganic carbon (calcium carbonate) using atomic absorption, or inductively-coupled plasma atomic absorption. What follows is a description of the methods for determining volume scattering, and ultimately backscattering, using a bench-top instrument, the Wyatt Technologies LASER light-scattering photometer. The methods are designed to estimate volume scattering of total particulate matter as well as particulate inorganic carbon ( $\text{CaCO}_3$  which involves measuring the VSF of seawater samples before and after sample acidification). These VSF measurements can be made in discrete water samples (Sect. 4.3), and along ship tracklines in a flow-through setup (Sect. 4.4).

---

<sup>3</sup> Certain commercial equipment, instrument, or materials are identified in this document to foster understanding. Such identification does not imply recommendation or endorsement by the National Aeronautics and Space Administration, nor does it imply that the materials or equipment identified are necessarily the best available for the purpose.

## 4.2 VOLUME SCATTERING FUNCTION INSTRUMENTS AND CALIBRATION

A Wyatt Technologies Dawn-F LASER Light-Scattering Photometer is used for measurement of the volume scattering function at 15 angles. It has a vertically polarized LASER source, and the particular instrument used for the work reported here has been fitted, at different times, with a Helium-Neon LASER (632.8 nm) and an Argon Ion LASER (514.0 nm). We have also used the newer Wyatt instrument, the EOS LASER light scattering photometer, equipped with a frequency-doubled Neodinium Yttrium Vanadate LASER (532 nm). The EOS instrument has a similar optical lay-out to the Dawn-F, and here we describe only the Dawn F instrument.

Figure 4.1 (Wyatt 1992) shows the physical layout of the LASER light source and detectors. The vertically-polarized LASER beam travels through a “doughnut” shaped, anodized aluminum block, with radial holes bored to the center “well”, which holds the glass cuvette. The instrument is equipped with two photodiodes for LASER power monitoring (one prior to passage through the viewing cuvette, and one post-passage). The detector ports are at the outside of the “doughnut”, and volume scattering is measured in a plane containing the LASER axis. For discrete measurements, the cuvette is cylindrical, and is inserted into the center of the well. For discrete aqueous measurements, the instrument viewing angles are 26.56°, 29.05°, 32.00°, 35.54°, 39.80°, 45.00°, 51.34°, 59.04°, 68.20°, 78.69°, 90.00°, 101.31°, 111.80°, 120.46°, 128.66°, 135.00°, 140.19°, and 144.46°. These angles are chosen to provide relatively small scattering angles for various liquid solvents, and to space the cotangents of the measurement angles at equal intervals. The first three detectors are useable for some solvents, but owing to the viewing geometries, and refraction by water and glass acting together with cell geometry, they cannot be used for measurements of aqueous samples. Complete scans of the VSF at the above angles are measured at 200 Hz, and averaged over 10 s. Examples of batch-mode data measured with this instrument can be found in Balch *et al.* (2001).

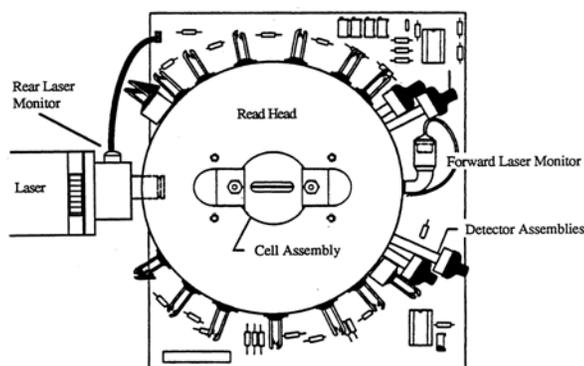


Figure. 4.1: Top view of the read head of the Dawn-F LASER light scattering photometer, showing the LASER and multiple detectors arranged radially around the central viewing volume. [Figure taken from the instrument manual (Wyatt 1992)].

The LASER beam in the Dawn-F has a  $e^{-2}$  Gaussian beam profile radius of 0.39 mm, which makes the effective viewing volume of the light-scattering photometer  $\sim 0.25$  mL, although this varies for each detector since the scattering volume is the intersection of the beam cross section and each detector’s field of view. The LASER beam divergence is 2.04 mrad, its polarization ratio is  $>500:1$  and optical noise is  $<1\%$ . Power stability of the Dawn is  $\pm 5\%$ . In terms of static alignment, the concentricity is  $<0.05$  mm and parallelism is  $<0.6$  mrad. For making discrete scattering measurements, a glass cuvette is placed in the center well of the instrument (Fig. 4.2). Glass scintillation vials are of the correct size to fit within the instrument, but extra care must be taken if they are used due to irregularities in the thickness of the glass walls. In a batch of 100 vials, for example, only 2 or 3 may have walls with minimal irregularities and thickness variations that are optically suitable for their use in these measurements.

Even when such vials are carefully selected for optical quality, they should be rotated slightly between replicate measurements to average the effects of surface irregularities. The best solution is to custom fabricate cuvettes from optical-grade glass, for use in this instrument.

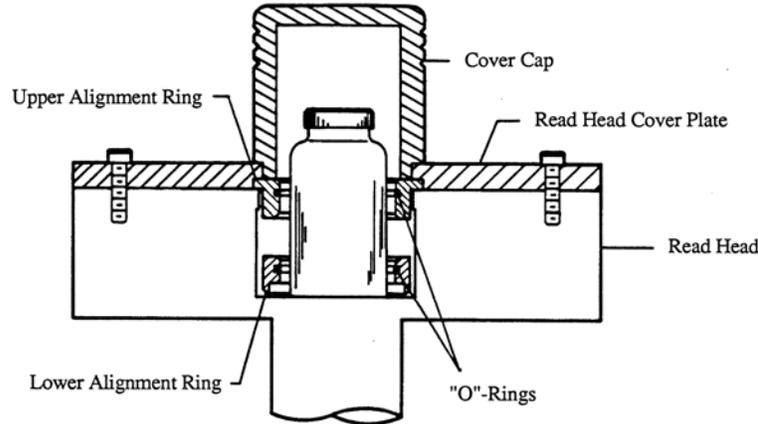


Figure 4.2: Side view of Dawn-F light scattering photometer read head, configured with a glass vial to make measurements of volume scattering in discrete water samples. [Figure taken from the instrument manual (Wyatt 1992)].

The Wyatt Technologies Dawn F Laser Light-Scattering Photometer can also be used for flow-through measurements. Most specifications are the same as given above for batch mode, except for the cuvette, which is adapted from one used for high-performance liquid chromatography applications. The flow-through cuvette is shown schematically in Figure 4.3. In this case, water flows through a polished glass cuvette, along the axis of the LASER beam, away from the LASER. In this manner, the cuvette walls are not near to the detected volume, which minimizes interaction of the LASER beam with cuvette walls. The “cuvette” is actually a cylindrical cavity in a glass disk, flattened at each end of the flow tube, that fills the read head enclosure so that the detectors view its glass-air interface at normal incidence; this geometry eliminates a second refractive direction change of the measured radiant intensities. Flow enters and exits the sensing volume through fittings outside the plane of VSF measurements. By virtue that the refractive index of the water is less than the cuvette glass, it is possible to measure the VSF at relatively small angles with low background noise. For flow-through measurements, the angles are slightly different than they are in batch mode, due to the composition and geometry of the cuvette. The 15 VSF angles ( $\psi$ ) are: 21.7°, 28.7°, 36.1°, 44.5°, 54.0°, 64.9°, 77.1°, 90.0°, 102.9°, 115.1°, 126.0°, 135.5°, 143.9°, 151.3°, and 158.3°.

The angular dependence of scattered light intensity measured using the Dawn-F instrument is expressed in terms of the **Rayleigh ratio**, defined as:

$$R(\psi) = \frac{I(\psi)r^2}{I(0)V}, \quad (4.1)$$

where  $I(\psi)$  is the scattered radiant intensity,  $I(0)$  is the radiant intensity of the incident beam,  $V$  is the volume of the intersection of the LASER beam and detector field-of-view, and  $r$  is the distance from the scattering volume to the detector. To determine the combined Rayleigh ratio for suspended particles  $R_p(\psi)$  and dissolved materials  $R_d(\psi)$  in water (or another liquid), the baseline scattered radiant intensity, e.g.  $I_w(\psi)$  due to pure water, may be subtracted to yield

$$R_p(\psi) + R_d(\psi) = \frac{[I(\psi) - I_w(\psi)]r^2}{I(0)V}, \quad (4.2)$$

where  $I_w(\psi)$  is determined by measuring the scattered radiant intensity distribution of optically pure water (Volume IV, Chapter 3) and normalizing it to remove differences between  $I(0)$  during the two measurements. The angular distribution of scattering by particles could be similarly separated from that associated with dissolved material by filtering the water samples through a  $0.2 \mu\text{m}$  filter, measuring the scattered radiant intensity distribution of the filtrate, and subtracting both  $I_w(\psi)$  and  $I_d(\psi)$  in an equation similar to (4.2); for seawater samples. In the present context, however, the scattering contribution of dissolved substances is typically negligible in all but river-dominated waters.

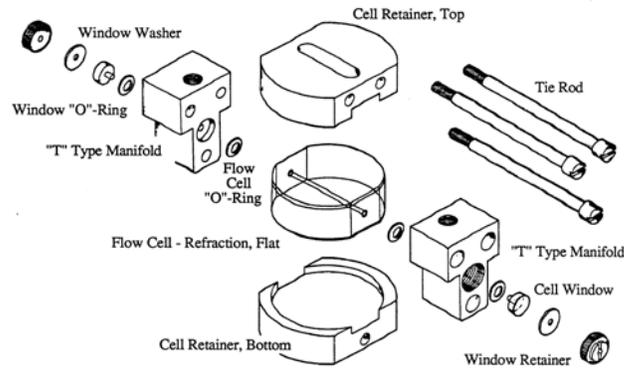


Figure 4.3: Exploded view of the Dawn F flow-through cuvette. Laser light passes along the axis of the flow cell. [Figure taken from the instrument manual (Wyatt 1992)].

The Wyatt Technologies Dawn Laser Light-Scattering Photometer is absolutely calibrated at  $\psi = 90^\circ$  using a solid isotropic scattering standard (the composition of the standard is proprietary to Wyatt Technologies; only calibration coefficients are available from the manufacturer). The solid standard gives a higher scattering signal than most organic solvents, also used in calibration of flow-through measurements (see below). As long as the solid standard is cleaned well on the exterior, it provides excellent consistency between calibrations, and the central sensed volume is always “dust free”. The Rayleigh ratio of the particular glass standard used by the authors, for example, is  $1.22 \times 10^{-4} \text{ cm}^{-1}$ .

Dark offsets are recorded during each calibration and used to correct measurements as

$$R(90^\circ) = A_{\text{CS}} \frac{S_{90} - S_{90}^{\text{dark}}}{S_0 - S_0^{\text{dark}}}, \quad (4.3)$$

where  $A_{\text{CS}}$  is the *configuration-specific* calibration constant (provided by the manufacturer),  $S_{90}$  and  $S_{90}^{\text{dark}}$  are the signal responses of the  $90^\circ$  detector with the LASER on and off, respectively, while  $S_0 - S_0^{\text{dark}}$  are the voltages of the LASER monitor detector, again with LASER on and off, respectively. To correct for specific instrument differences and geometrical factors that may be unique to each instrument, a true instrument calibration constant,  $A_{\text{inst}}$ , is defined as

$$A_{\text{inst}} = A_{\text{CS}} \frac{F}{n_L n_G}, \quad (4.4)$$

where  $n_L$  and  $n_G$  are the liquid and cuvette refractive indices, respectively, and  $F$  represents Fresnel reflection losses at the various interfaces in the cell as

$$F = \left[ 1 - \left( \frac{n_G - n_L}{n_G + n_L} \right)^2 \right]^2 \left[ 1 - \left( \frac{n_G - 1}{n_G + 1} \right)^2 \right]. \quad (4.5)$$

With the two calibration constants,  $A_{inst}$  and  $A_{CS}$ , the Dawn  $F$  can be calibrated with one solvent, and the results of that calibration may be applied to measurements made with another solvent.

Detector signals are normalized to the calibrated Rayleigh ratio at  $90^\circ$  using the above-mentioned isotropic glass scattering standard, to correct for the fact that each detector views a different scattering volume and subtends a different solid angle with respect to the scattering volume.

Characterization of the Wyatt in flow-through mode is identical to batch mode, except for the cuvette (see above). In flow-through mode, normalization of the detectors is achieved by running a solution of Dextran (Mol. Wt 39,200; Sigma 9004-54-0) through the flow-through cuvette. This solution scatters isotropically (Wyatt 1992). Absolute calibration of the instrument is achieved by pumping  $0.02\mu\text{m}$  filtered, analytical grade methanol, toluene, or optically pure water through the optical cell. The software can accommodate particle-free methanol or toluene as a calibration standard. Toluene has the highest Rayleigh ratio ( $1.406 \times 10^{-5} \text{ cm}^{-1}$  at  $632.8\text{nm}$ ) of any of the commonly-available solvents, and is therefore useful for calibration. The Rayleigh ratios for various solvents and wavelengths are embedded in the instrument software.

Quality control of the Wyatt Technologies measurements is checked by viewing  $0.02\mu\text{m}$ -filtered, Milli-Q water and comparing the VSF at  $90^\circ$  and derived  $b_b$  to published values for pure water (Vol. I, Ch. 2 and Vol. IV, Ch. 1). For ultra filtered, HPLC-grade distilled water measurements, our software fits a standard Rayleigh function to the volume scattering data instead of the Beardsley-Zaneveld (1969) polynomial function (which is fitted to volume scattering measurements of suspended particles). Accurate calibration is evident immediately when viewing the fitted volume scattering function along with actual data.

### 4.3 DISCRETE WATER SAMPLES: SAMPLE ACQUISITION, TREATMENT, AND VSF MEASUREMENTS

#### *Sample Acquisition*

Discrete seawater samples for VSF measurements can be drawn from a bucket, or from Niskin bottles, provided that samples are well-mixed, with particles suspended. It should be noted that if Niskin bottles are left on deck without stirring, large, negatively buoyant, mineral particles can settle below the sampling valve. The approximate volume needed for discrete measurements with the Dawn  $F$  is 10 mL to 15 mL. Prior to running discrete seawater samples, vials are checked for any irregularities in wall thickness, and blanks are run using  $0.02 \mu\text{m}$ -filtered, HPLC-grade distilled water. In the event that particles do not remain in suspension over the period of a measurement, samples can be regularly swirled, or a small magnetic stirring motor can be installed below the sample well and a thoroughly-cleaned magnetic stir bar (“flea”) placed within the sample vial. Care should be taken to slightly rotate the sample cuvette between sample replicates to average the effects of any irregularities in the walls of the cuvette.

#### *VSF Measurements of Untreated Seawater Samples*

Volume scattering measurements of discrete, or “batch”, samples are performed using 3 to 5 replicates of 10 s measurements sampled at 200 Hz; this yields to 6000-10000 VSF measurements, which are averaged. The replicated measurement sets are combined to determine the average total backscattering coefficient  $b_b(\lambda)$  (Sect. 4.6 below) for the sample.

#### *Sample Acidification to Dissolve Particulate $\text{CaCO}_3$*

To acidify discrete samples and dissolve  $\text{CaCO}_3$  coccoliths, seawater aliquots can be either bubbled with  $\text{CO}_2$  for 30 s, or  $0.02 \mu\text{m}$ -filtered acetic acid may be added (0.05 % final concentration) to give a final pH < 5.8. Acid addition is preferred over  $\text{CO}_2$  bubbling, because the bubbling process can induce formation of more particles.

*VSF Measurements of Acidified Samples*

Following acidification to dissolve particulate  $\text{CaCO}_3$ , 3 to 5 additional replicate 10 s VSF measurements are recorded at 200 Hz. The overall set of VSF measurements is then combined to determine the average “acidified” backscattering coefficient  $b_b^{\text{acid}}(\lambda)$  of the seawater sample.

*Quality Control and Uncertainty*

Quality control is assured by frequent checks of the volume scattering of 0.02 $\mu\text{m}$  filtered, Milli-Q water. Based on more than 5 years of measurements of “particle-free” water with the Wyatt instrument, the Type A uncertainty of the VSF determinations at 632.8 nm is approximately  $1.5 \times 10^{-4} \text{ m}^{-1} \text{ sr}^{-1}$ .

## 4.4 CONTINUOUS FLOW-THROUGH MEASUREMENTS ALONG SHIP TRACKS: SAMPLE ACQUISITION, TREATMENT, AND VSF MEASUREMENTS

*Flow-Through Sampling Apparatus and Debubbler*

Water flow in the continuous underway system is first run through three de-bubblers, arranged in series, to remove small bubbles (and their associated light scattering). A metering pump then delivers the water to the Wyatt Technologies Dawn F light scattering photometer at an approximate flow rate of 11  $\text{mL min}^{-1}$ . It is easy to detect bubble contamination either by visual inspection of the cuvette during measurements, or retrospectively, by screening the data for evidence of the high scattering levels that occur when bubbles are present in the LASER beam.

*In Line Acidification and Flushing Method*

To estimate the backscattering of particulate  $\text{CaCO}_3$  in raw seawater measurements, a peristaltic pump is activated, following VSF measurements of raw seawater, to inject 0.05 % (final concentration) glacial acetic acid into the flow stream. Flow through a mixing coil insures adequate mixing of the acid and seawater. A micro-pH probe downstream of the optical cuvette monitors the pH. The conventional solubility product for calcite in seawater at 25° C and 35 PSU is  $K_{\text{sp}}^{\text{calc}} = 4.467 \times 10^{-7}$  (UNESCO 1987). This value assumes that the activity of the pure solid phase is unity, and that the calcite is in its chemically pure form (Stumm and Morgan 1981); these are both reasonable assumptions. When the pH of the sea water flow is  $< 5.8$  (since  $\text{p}K_{\text{sp}}^{\text{calc}} = 6.35$ , a pH less than this value assures dissolution of  $\text{CaCO}_3$ ), the VSF of the acidified seawater flow is ready to be measured, without scattering contributions from particulate  $\text{CaCO}_3$ .

*Measurement Cycles: Untreated Sample VSF – Acidified Sample VSF – System Flushing*

In flow mode, the Wyatt Technologies Model Dawn LASER light scattering photometer has seawater flowing at an approximate rate of 11  $\text{mL min}^{-1}$ . As with the batch measurements, all detectors can be scanned at rates up to 400 Hz, but for most flow-through applications we slow the scanning rate to 200 Hz in order not to sample the same seawater volume twice.

Labview software is used to control all aspects of sampling. The VSF, at all angles specied above (Sect. 4.2) for the flow-through configuration, are measured at 200 Hz, and averaged over 1 s intervals for calculation of the backscattering coefficient. The 1 s averages are recorded for 50 s, and combined to determine the average total backscattering coefficient  $b_b(\lambda, t)$  (Sect. 4.6 below). This measurement cycle is first performed to determine the VSF, and  $b_b(\lambda, t)$  for raw seawater. Following acidification to dissolve particulate  $\text{CaCO}_3$ , data from the Wyatt Dawn light scattering photometer are recorded and averaged for 50 s to determine  $b_b^{\text{acid}}(\lambda, t)$ . Finally, the acid pump is stopped, and the flow-through system is flushed until the pH returns to the alkaline values of raw seawater. The entire measurement cycle is then repeated.

The time for a complete acidification cycle can be adjusted, but we have preferred to collect average VSF measurements such that one complete raw/acidification cycle takes 4 min. At a typical ship speed of 10 knots ( $5 \text{ m s}^{-1}$ ), a 4 min measurement cycle represents an alongtrack distance of 1200 m, which is comparable to the resolution of satellite ocean color measurements using, *e.g.*, SeaWiFS and MODIS.

### *Quality Control*

Quality control is assured by frequent checks of the measured VSF of  $0.02 \mu\text{m}$  filtered, Milli-Q water. The flow-through cuvette can be observed through a viewing port on top of the cuvette to check for bubbles or bio-fouling. Essentially, the sides of the flow tube become visible if any biofouling occurs. If this happens, then the cuvette should be rinsed, and if necessary, disassembled and cleaned.

Routinely, at intervals of 24 hours, the flow is stopped and the flow cell is cleaned to remove bio-fouling. Storage of the cell with ultra clean solvent, or detergent, solves most problems associated with wall coatings by organisms. Sonication of the cuvette also helps maintain cleanliness. Complete cleaning requires removal of the end windows, and cleaning of the internal bore with solvent soaked lens paper. To correct for any bio-fouling, or instrument drift,  $0.02 \mu\text{m}$  filtered Milli-Q distilled water is pushed through the flow cell to adjust the calibration at frequent intervals throughout the duration of a measurement sequence. After the cruise, all data are corrected using the distilled water calibration results, by comparison to the known backscattering of pure water.

LASER alignment, which is a critical characteristic for good instrument performance, and has proven to be extremely stable with the Dawn F.

As with the discrete measurements above, based on water measurements of “particle-free” water, measurements of the VSF with the Wyatt Dawn in flow mode have a Type A uncertainty of  $1.5 \times 10^{-4} \text{ m}^{-1} \text{ sr}^{-1}$ . This combined standard uncertainty includes uncertainties associated with instrument calibration and bio-fouling of the flow-through cuvette during continuous use.

## 4.5 ANCILLARY MEASUREMENTS

### *Coccolith and Coccolithophore Enumeration*

Microscope enumeration of coccolithophores and coccoliths is done by filtering a 50 mL water sample through a Millipore HA filter. The filter is then rinsed with borate buffer and frozen in a Petri dish until the time that the cells and coccoliths will be counted (Haidar *et al.* 2000; Haidar and Thierstein 2001). To prepare the sample for counting, the filter is placed on a glass microscope slide, and  $60^\circ\text{C}$  Canada Balsam is placed on top of the filter, followed by a cover slip. The clarified filter is examined with an Olympus BH2 microscope equipped with polarization optics. Coccoliths and plated coccolithophores can then be counted based on their unique birefringence patterns. For an example of these patterns, see Moshkovitz (1989). For statistical reasons, 200 coccoliths or cells are counted from each sample, when that number or more are present on the filter.

### *Particulate Inorganic Carbon*

Particulate inorganic carbon (PIC) is estimated from measurements of particulate calcium. Samples of 100 mL to 300 mL are filtered through  $0.4 \mu\text{m}$  poresize polycarbonate filters. The filters are then rinsed with filtered seawater and borate buffer ( $\text{pH} = 8$ ) to remove  $\text{CaCl}_2$  (Fernández *et al.* 1993), and are placed in centrifuge tubes with 5 mL of 0.5 % Optima grade  $\text{HNO}_3$ . Particulate Ca is measured according to Balch, Drapeau, and Fritz (2000), except that an inductively-coupled plasma atomic absorption spectrometer is now used to measure calcium, instead of a graphite furnace atomic absorption spectrometer. The sensitivity of the technique, after correction for the volume of seawater filtered, is about  $2 \text{ ng Ca L}^{-1}$ . The coefficient of variation for the measurements is  $\sim \pm 3 \%$  at a PIC concentration of  $\sim 1 \mu\text{gC L}^{-1}$ . By using polycarbonate filters, any Ca contamination from residual, interstitial water in the filter is minimized. This helps lower blank values.

## 4.6 DATA ANALYSIS METHODS

Backscattering is calculated by integrating the VSF measurements from 90° to 144.5° using standard trapezoidal integration. Integration of backscattering from 144.5° to 180° is accomplished by fitting a polynomial function (Beardsley and Zaneveld 1969) to VSF measurements at 45°, 90°, and 135°. The polynomial function is then integrated, and the two integrals (90° to 144.5° and 144.5° to 180°) are summed to estimate backscattering coefficients  $b_b(\lambda)$  and  $b_b^{\text{acid}}(\lambda)$  from the raw and acidified seawater samples, respectively. The difference between integration using the above method, versus simply integrating the fitted polynomial between 90° and 180°, shows differences < 5%, in accordance with the observations of Gordon (1976; his Table 1). Finally, the backscattering coefficient for particulate CaCO<sub>3</sub> (calcite) is calculated as

$$b_b^{\text{calc}}(\lambda) = b_b(\lambda) - b_b^{\text{acid}}(\lambda) \text{ m}^{-1}. \quad (4.6)$$

The values of calcite backscattering can be used to estimate the quantity of suspended calcium carbonate, provided one has an estimate of the backscattering cross-section of coccoliths (Balch *et al.* 1999). Discrete estimates of particulate inorganic carbon can be used to calibrate the optically-derived estimates.

## 4.7 DISCUSSION: PROTOCOL STATUS AND FUTURE DIRECTIONS

The first results of this instrument were published in 1999 (Balch *et al.* 1999), in which a flow cytometer was used to sort individual calcium carbonate particles into the cuvette. Continuous, surface  $b_b$  observations from the Arabian Sea have also been described (Balch *et al.* 2001).

The Wyatt Technology Dawn photometer and HOBILABS Hydroscat 2 scattering photometer are plumbed in series for our NASA Gulf of Maine ferry program (Fargion and McClain 2001). Both instruments are used to continuously log backscattering along a 325 Km track between Yarmouth, Nova Scotia and Portland, ME. This has provided a large data set from which to compare the two instruments. Particulate backscattering measured with the Hydroscat-2, has been compared to particulate backscattering measured with the Wyatt Technology instrument. To do these comparisons our Hydroscat-2 was directed into a 21 L volume of water and  $b_b$  estimates were based on a VSF measurement at one angle and two wavelengths (assuming one shape for the VSF). The Wyatt instrument estimated average  $b_{b \text{ pg}}$  in 9 mL to 12.5 mL of seawater by measuring the VSF at 15 angles and one wavelength at 200Hz (such  $b_b$  calculations allowed the shape of the VSF to vary). Data from our 1999 and 2000 field seasons were used for this comparison. The Hydroscat 2 data at 476 nm and 676 nm were interpolated to estimate the 514 nm value. The interpolation assumed that backscattering increases as the wavelength raised to a power (Mobley 1994). The results (Fig. 4.4) produced a cloud of data approximately centered about the 1:1 line, but with significant variability. Generally, the two instruments tracked each other quite well, with divergences occurring primarily in very turbid water. The best-fit relationship between the two instruments was:  $b_{\text{bp}514 \text{ Hobi}} = 0.19 (b_{\text{bp}514 \text{ Wyatt}})^{0.715}$  [ $r^2 = 0.423$ ;  $n = 3029$ ;  $P < 0.001$ ]. Some of the differences can be ascribed to the shape of the VSF implicitly assumed in the HOBILABS instrument. Detailed comparisons between VSF at the same angles, as measured by the HOBILABS and Wyatt Technology instruments, has been published elsewhere (Vaillancourt *et al.* 2003).

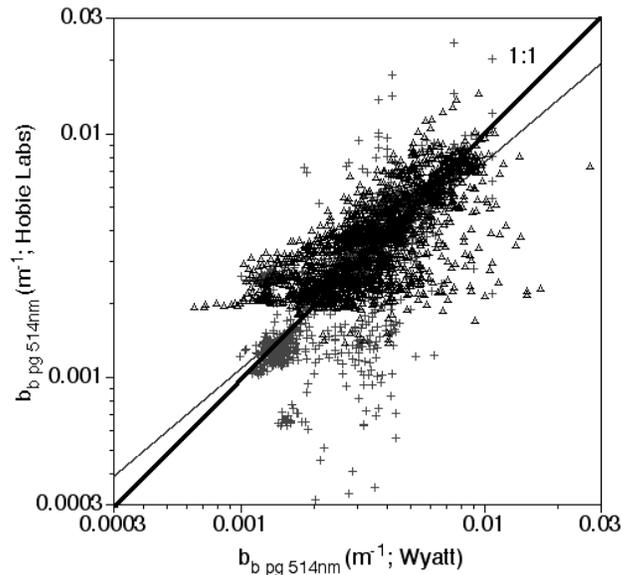


Fig. 4.4: Comparison between particulate backscattering estimates by the Wyatt Dawn F LASER light scattering photometer and the HOBILABS Hydroscat-2 (see text). Data from 1999 (+ symbols) and 2000 (triangles), continuous underway sampling in the Gulf of Maine were used in this comparison.

## ACKNOWLEDGEMENTS

Special thanks go to Bruce Bowler (Bigelow Laboratory for Ocean Sciences) and Charlie Byrne (Univ. Miami), who wrote our software for collecting data from the Wyatt Dawn F and EOS light scattering photometers while at sea. Thanks also go to Katherine Kilpatrick (Univ. Miami) for her earlier help in setting up the Dawn F scattering photometer for shipboard use. Dr. Jennifer Fritz (Univ. Miami), Amanda Ashe (Oregon State University) and Emily Booth (Bigelow Laboratory for Ocean Sciences) helped collect shipboard data sets on volume scattering. Dr. Ken Voss (Univ. Miami) helped us perform an intercalibration of various light scattering photometers, including the Wyatt instruments described here. Lastly, we'd like to thank the many individuals at Wyatt Technology Corporation (Santa Barbara, CA, U.S.A.) for their helpful advice on using their scattering photometers for oceanographic applications. Methods development, instrumentation, and observations described herein were generously supported by various grants to W.M.B. over the years: National Aeronautics and Space Administration (NAGW2426; NAS5-97268; NAS5-31363; NAG5-10622; NASA EPSCOR EP-02-14); Office of Naval Research (ONR N00014-91-J-1048, N00014-97-1-0034; N00014-98-1-0882; N00014-99-1-0645; N00014-01-1-0042), National Science Foundation (OCE-9022227; OCE-9596167), and NOAA (NA56RM0258; 40-AA-NE-005996).

## REFERENCES

- Balch W.M., D. Drapeau, and J. Fritz, 2000: Monsoonal forcing of calcification in the Arabian Sea. *Deep-Sea Research II*, **47**: 1301-1337
- Balch W.M., Drapeau D., J Fritz, B. Bowler, and J. Nolan, 2001: Optical backscattering in the Arabian Sea-continuous underway measurements of particulate inorganic and organic carbon. *Deep Sea Research I*, **48**: 2423-2452

- Balch W.M., D.T. Drapeau, T.L. Cucci, R.D. Vaillancourt, K.A. Kilpatrick, and J. Fritz, 1999: Optical backscattering by calcifying algae--Separating the contribution by particulate inorganic and organic carbon fractions. *J. Geophys. Res.* **104**: 1541-1558
- Balch W.M., R.W. Eppley, M.R. Abbott, and F.M.H. Reid, 1989: Bias in satellite-derived pigment measurements due to coccolithophores and dinoflagellates. *J. Plankton Res.*, **11**: 575-581
- Beardsley, G.F. and J.R.V. Zaneveld (1969) Theoretical dependence of the near-asymptotic apparent optical properties on the inherent optical properties of sea water. *J. Opt. Soc. Am.*, **59**: 373-377
- Fargion, G.S. and C.R. McClain, 2001: *SIMBIOS project 2000 annual report*. National Aeronautics and Space Administration, Goddard Space Flight Center, Greenbelt, MD
- Fernández, E., P. Boyd, P.M. Holligan and D.S. Harbour, 1993: Production of organic and inorganic carbon within a large scale coccolithophore bloom in the northeast Atlantic Ocean., *Mar. Ecol. Prog. Ser.*, **97**: 271-285
- Gordon, H.R., 1976: Radiative transfer in the ocean: A method for determination of absorption and scattering properties., *Appl. Opt.*, **15**: 2611-2613
- Haidar, A.T. and H.R. Thierstein, 2001: Coccolithophore dynamics off Bermuda (N. Atlantic). *Deep Sea Research*, **48**: 1925-1956
- Haidar, A.T., H.R. Thierstein and W.G. Deuser, 2000: Calcareous phytoplankton standing stocks, fluxes and accumulation in Holocene sediments off Bermuda (N. Atlantic). *Deep Sea Research*, **47**: 1907-1938
- Mobley, C.D. 1994: *Light and water: Radiative transfer in natural waters*. Academic Press, New York
- Moshkovitz, S. and K. Osmond, 1989: The optical properties and microcrystallography of Arkhangelskiellaceae and some other calcareous nannofossils in the Late Cretaceous. In: Crux, J.A. and S.E. van Heck (eds) *Nannofossils and their applications*. Ellis Horwood, Chichester, pp 76-97
- Stumm, W., J.J. Morgan, 1981: *Aquatic Chemistry*. John Wiley & Sons, Inc., New York
- UNESCO (1987) *Thermodynamics of the carbon dioxide system in seawater. Report by the carbon dioxide sub-panel of the joint panel on oceanographic tables and standards*. United Nations Educational, Scientific and Cultural Organization, 51, Paris
- Vaillancourt, R.D., C.W. Brown, R.R.L. Guillard and W.M. Balch, 2003: Light backscattering by phytoplankton algae and a heterotrophic bacterium measured with a multi-wavelength, fixed-angle backscattering meter. *Journal of Plankton Research* (in press)
- Wyatt, P., 1992: *Instruction manual for the DAWN Model F.*, Wyatt Technology Corporation, Santa Barbara, CA